

August 2011

Single-Analyte ELISArray for Rat Handbook

For quantitative measurement of the amount of an individual cytokine or chemokine using ELISA in multiple samples



Sample & Assay Technologies

QIAGEN Sample and Assay Technologies

QIAGEN is the leading provider of innovative sample and assay technologies, enabling the isolation and detection of contents of any biological sample. Our advanced, high-quality products and services ensure success from sample to result.

QIAGEN sets standards in:

- Purification of DNA, RNA, and proteins
- Nucleic acid and protein assays
- microRNA research and RNAi
- Automation of sample and assay technologies

Our mission is to enable you to achieve outstanding success and breakthroughs. For more information, visit www.qiagen.com.

Product Use Limitations

Single-Analyte ELISArray Kits for Rat are intended for molecular biology applications. These products are not intended for the diagnosis, prevention, or treatment of a disease.

CONTENTS

I.	Background and Introduction	4
II.	Materials Provided	6
III.	Precautions	6
IV.	Additional Materials Required	6
V.	Complementary Products	6
VI.	Protocol	7
	A. Sample Preparation and Handling	7
	B. Reagent Preparation	8
	C. Assay Procedure	9
	D. Data Analysis	11
VII.	Troubleshooting and Frequently Asked Questions	11

I. Background and Introduction

The-Single Analyte ELISArray Kits are designed to quantitatively measure the amount of an individual protein analyte using a standard sandwich enzyme-linked immunosorbant assay (ELISA) technique. A target-specific capture antibody has been coated onto 8-well strips placed into a 96-well plate format. In addition to the pre-coated plate, the kit also includes the detection antibody, Antigen Standard, and a complete set of reagents for a colorimetric ELISA. The performance of other ELISArray Kits currently on the market is limited by the quality of the antibody pair produced by their own company. QIAGEN's ELISArray Kits are developed with the best possible antibody pairs on the market. To develop these high-performance ELISArray Kits, several antibody pairs were screened by comparing them side-by-side on signal intensity, specificity, and antigen response dynamic range. The best possible pair was chosen for each ELISArray Kit. Take advantage of SABiosciences' assay development and optimization expertise with the Single-Analyte ELISArray Kits.

The Single-Analyte ELISArray Kits are quantitative sandwich-based enzyme-linked immunosorbant assays. Figure 1 displays an overview of the assay protocol. A capture antibody specific for a single protein has been coated onto the 8-well strips. Replicate serial dilutions of a standard for the protein of interest are added to one set of replicate strips, while replicate dilutions of unknown biological samples are added to other sets of strips. Incubation allows the capture antibody to bind the specific protein of interest. After washing away unbound protein, biotinylated detection antibody added to the wells also binds the captured analyte. After washing again to remove unbound material, an avidin-horseradish peroxidase conjugate is added. The wells are again washed and a colorimetric substrate solution is added, which produces a blue color in direct proportion to the amount of protein analyte present in the initial sample. The color development is stopped by adding stop solution, and the absorbance at 450 nm is read. The readings from the unknown samples are then compared to the standard curve to quantify the amount of protein in the original sample.

Benefits of the Single-Analyte ELISArray Kits:

- **Verified High-Performance Assays:** Low background. High sensitivity. Excellent linearity.
- **Easy to Use:** No special equipment required, only the ELISA plate reader already in your lab.

Figure 1: Overview of the Single-Analyte ELISA Kit Procedure.
Total Time ~ 4.5 hours

1. Prepare all reagents.
Set up experimental samples, Antigen Standard, and negative control.



2. Add 50 μ l assay buffer into each well of ELISArray plate.
Transfer 50 μ l of samples and serial diluted Antigen Standard into the appropriate wells of the ELISArray plate.
Incubate 2 hours.



3. Wash three times.



4. Add 100 μ l Detection Antibody Solution. Incubate 1 hour.



5. Wash three times.



6. Add 100 μ l Avidin – HRP. Incubate 30 minutes.



7. Wash four times.



8. Add 100 μ l Development Solution. Incubate 15 minutes in the dark.



9. Add 100 μ l Stop Solution. Read OD 450 nm within 30 minutes.

II. Materials Provided:

Component / Description	Quantity
BOX 1	
Antigen Standard (1 µg/ml)	1.5 ml tube
Detection Antibody	1.5 ml tube
Avidin-HRP Conjugate	1.5 ml tube
10% BSA	15 ml bottle
BOX 2	
Capture Antibody-Coated 8-well strips	One plate with 12 x 8-well strips
Sample Dilution Buffer Stock	60 ml bottle
Assay Buffer Stock	60 ml bottle
Wash Buffer (10x Concentrate)	125 ml bottle
Development Solution	25 ml bottle
Stop Solution	60 ml bottle

Storage Conditions:

Box 1 is shipped on dry ice or blue ice packs and should be stored in **non-frost-free freezer** at -20°C upon receipt.

Box 2 is shipped at ambient temperature and should be stored at 4°C upon receipt.

Do not use kit beyond the expiration date printed on the label.

Store any unused 8-well strips from an opened pouch in a zip-lock bag at 4°C.

III. Precautions

1. The Development Solution is toxic if inhaled or swallowed. Avoid contact with skin. Keep container tightly closed when not in use.
2. Stop Solution is an acidic solution. Wear eye, hand, face, and clothing protection when using this material.

IV. Additional Materials Required:

1. Standard ELISA Microplate Reader
Capable of measuring 450 nm absorbance with a 570 nm correction wavelength
2. Calibrated Multi-Channel Pipettor
3. Wash Bottle
4. Microcentrifuge Tubes
5. Laboratory Timer

V. Complementary Products:

Multi-Analyte ELISArray Kits for Rat

VI. Protocol:

Please read through this entire protocol before beginning your experiment.

A. Sample Preparation and Handling:

High quality input material is **ESSENTIAL** for obtaining good results.

The most important prerequisite for any ELISA analysis experiment is consistent, high-quality experimental samples. Therefore, the sample preparation and handling procedures are critical to the success of the experiment. Specimens should be clear, because residual traces of particulate matter, heme, lipids or other contaminants will interfere with the performance of the ELISA.

Recommended Preparation Methods:

High quality protein samples for the ELISA experiment must be prepared using one of the following methods, each specific for your biological sample:

a. Cell Culture Supernatants:

Remove any particulate material by centrifugation for 10 minutes at 1000 x *g* and assay immediately, or aliquot and store samples at $\leq -20^{\circ}\text{C}$. Avoid repeated freeze / thaw cycles.

b. Serum:

Use a serum separator tube and allow samples to clot for 30 minutes before centrifugation for 15 minutes at 1000 x *g*. Remove serum and assay immediately or store samples at $\leq -20^{\circ}\text{C}$. Avoid repeated freeze / thaw cycles.

c. Plasma:

Collect plasma using EDTA as an anticoagulant. Within 30 minutes of collection, centrifuge for 10 minutes at 1000 x *g*. Assay immediately or store samples at $\leq -20^{\circ}\text{C}$. Avoid repeated freeze / thaw cycles.

Note: Rat serum samples require a 2-fold dilution with sample dilution buffer before proceeding with the assay.

B. Reagent Preparation:

Use only polypropylene tubes for the antigen and detection antibody steps.

Briefly centrifuge any 1.5 ml tube in the kit before opening it in order to collect its contents at the bottom of the tube.

1. ELISA Array Kit Reagents:

- a. Thaw the 10% BSA at room temperature. Store on ice.
- b. Bring the Wash Buffer Concentrate, Assay Buffer Stock, Sample Dilution Buffer Stock and ELISArray plate to room temperature. Do not remove the ELISArray plate from its pouch at this time.
- c. Once finished with each reagent, return it to its proper storage conditions.
- d. Keep all other kit components at their recommended storage temperature until recommended by the protocol.

2. Wash Buffer

Visually inspect the Wash Buffer Concentrate to ensure that all components are in solution. If any precipitates are visible, briefly shake the bottle to suspend any precipitate. Dilute 50 ml of Wash Buffer Concentrate into de-ionized or distilled water (dH₂O) to a final volume of 500 ml. Transfer to wash bottle. Keep at room temperature.

3. Assay Buffer

Dilute 0.6 ml of 10% BSA into a final volume of 30 ml with Assay Buffer Stock. Keep at room temperature.

4. Sample Dilution Buffer:

Dilute 2 ml of 10% BSA into a final volume of 20 ml with Sample Dilution Buffer Stock. Store on ice.

C. Assay Procedure:

NOTE: *Very accurate and precise pipetting is critical to the reliability and consistency of any ELISA experiment. Make sure that all of your single- and multi-channel pipettors are calibrated before starting this protocol. Maintain incubation times and temperatures as consistent as possible across strips and experiments for best results.*

1. Prepare Antigen Standard dilutions and sample dilutions:
 - a. Just before use, thaw the Antigen Standards on ice for 20 minutes. Vortex gently to mix well.
 - b. Prepare a serial dilution of the Antigen Standard (2000, 1000, 500, 250, 125, 61.5, and 31.25 pg/ml) in a set of eight 5 ml (12 x 75) polypropylene tubes as follows:
 - i. Add 2495 μ l of Sample Dilution Buffer to the first tube, and 500 μ l of the same buffer to each of the remaining seven tubes.
 - ii. Add 5 μ l of the Antigen Standard to the first tube to make 2000 pg/ml Antigen Standard. Mix well.
 - iii. Generate a series of 1:2 dilutions by transferring 500 μ l from the first to the second tube, mixing well and then repeating from the second to the third and so forth on up to the seventh tube, as demonstrated in Figure 2.
 - iv. Do not add any source of Antigen Standard to the eighth tube to setup a required assay performed in the absence of antigen to determine the background signal.

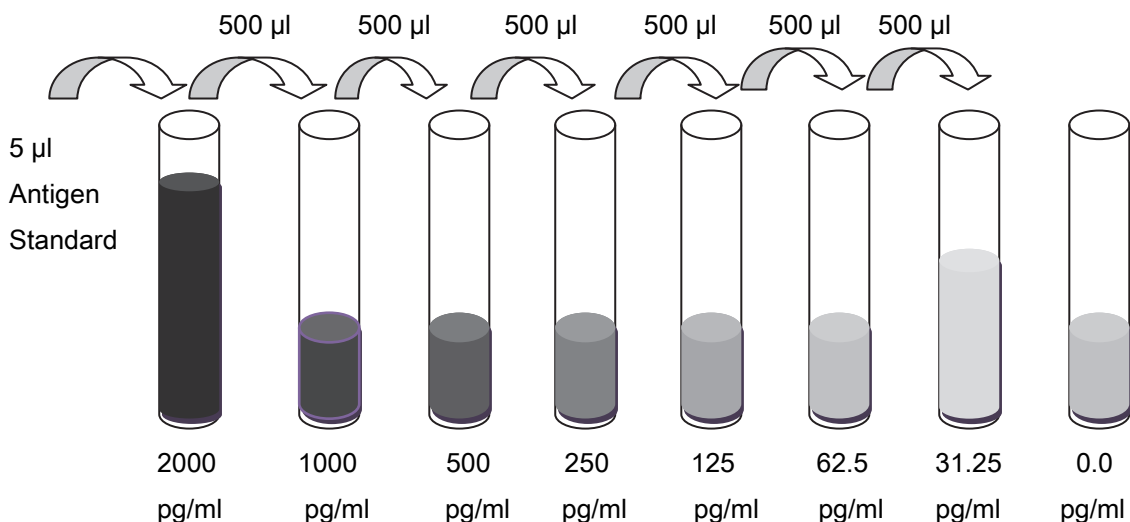


Figure 2: Generating a 1:2 Dilution Series and Negative Control for a Standard Curve.

- c. A series of dilutions of sample should be prepared in similar fashion using the same sample dilution buffer. Choose your serial dilution series based on the expected concentration of the analyte in your samples.

2. Remove the ELISArray plate from its pouch.
Using a multi-channel pipettor, add 50 μ l of Assay Buffer into each well of the ELISArray plate.

NOTE: *For best results, add all samples and reagents carefully by touching the pipette tips to the side of the wells just above the bottom and allowing the dispensed volume to run down the well wall. Do not blow out the last drop of volume dispensed to avoid introducing bubbles into the wells.*
3. Transfer 50 μ l of either the Antigen Standard or your sample serial dilutions to the appropriate wells of the ELISArray 8-well strips. Use multiple wells or strips to obtain replicate measurements.

NOTE: *Change your pipette tips with every addition to avoid cross-contamination of the wells.*
4. Cover the plate and gently shake or tap the strips for 10 seconds to mix.
Incubate for 2 hours at room temperature on your bench-top.
5. Warm the Development Solution and Stop Solution to room temperature in lab bench drawer protected from light.

NOTE: *Avoid prolonged exposure of the Development Solution to light or contact with water, air, or extreme temperature.*
6. Dilute the Detection Antibodies just before the next step:
Thaw the Detection Antibodies on ice for 30 minutes.
Dilute 0.55 ml of Detection Antibody to a final volume of 11 ml with Assay Buffer.
Mix well but gently.
7. Washing ELISArray Wells
Decant or aspirate the contents of the wells. Wash wells by filling with 350 μ l of 1x Wash Buffer. Gently shake or tap the plate for 10 seconds to mix and follow by decanting or aspirating the solution. Blot the strips upside down on absorbent paper to remove any residual buffer. Repeat twice for a total of three washes.
8. Add 100 μ l of Detection Antibody solution to each well.
Cover and gently shake or tap the plate for 10 seconds to mix.
Incubate 1 hour at room temperature on your bench-top.
9. Prepare Avidin-HRP:
Just before use, thaw the Avidin-HRP Conjugate on ice for 20 minutes.
Vortex gently to mix well.
Add 11 μ l of the Avidin-HRP Conjugate to 11 ml of Assay Buffer.

NOTE: *Avoid prolonged exposure of the Avidin-HRP Solution to light.*
10. Wash ELISArray wells as in step 7.
11. Add 100 μ l of Avidin-HRP to all wells.
Gently shake or tap the plate for 10 seconds to mix.
Incubate for 30 minutes at room temperature in the dark.
12. Wash ELISArray wells as in step 7, except for a total of 4 washes.

13. Just before use, transfer the needed amount of the Development Solution (100 μ l per well plus extra for viscosity pipetting errors) from its original bottle into a clean multi-channel pipettor reservoir.

NOTE: Use a different clean multi-channel reservoir for each solution to be dispensed.

14. Add 100 μ l of the Development Solution to each well.
Incubate the plate for 15 minutes at room temperature in the dark.
15. Add 100 μ l of Stop Solution to each well.
The color in the wells should change from blue to yellow.
16. Read absorbance at 450 nm within 30 minutes of stopping the reaction. If wavelength correction is available, subtract readings at 570 nm from the reading at 450 nm.

NOTE: Subtracting the A450 reading by the A570 reading corrects for any minor optical imperfections in the ELISArray plate. Uncorrected A450 readings may yield artificially high signals.

D. Data Analysis:

1. Processing the Raw Data

Typical absorbance values should range from 0.00 to 2.50. Calculate the mean (average) absorbance for each set of replicate standards and experimental samples. Subtract these values by the mean (average) background absorbance as determined from assays performed in the absence of antigen.

2. Standard Curve

- a. Plot the corrected mean absorbance for the standard on the y-axis against its concentration on the x-axis in a log-log graph for a seven-point standard curve.
- b. Calculate the best straight-line curve fit through the points on the graph within the linear dynamic range of the results. Do not include points at the low and high concentration extremes where the absorbance does not change greatly with differing concentration.

3. Determination of Unknowns

- a. Use the background-corrected mean absorbance of each dilution of each unknown experimental sample and the standard curve to determine the protein amount in the sample.
- b. If the experimental unknown samples have been diluted, the concentration read from the standard curve must be multiplied by the dilution factor.

VII. Troubleshooting and FAQs

For technical assistance and more information, please see our Technical Support Center at www.qiagen.com/Support or call one of the QIAGEN Technical Service Departments or local distributors (see back cover or visit www.qiagen.com).

Ordering Information

Product	Contents	Cat. no.
Single-Analyte ELISArray Kits for Rat	Antigen Standard, Detection Antibodies, Avidin-HRP Conjugate, 10% BSA, Donkey Serum, Precoated Capture Antibody 8-Well Strips, Sample Dilution Buffer Stock, Assay Buffer Stock, 10x Wash Buffer, Development Solution, Stop Solution	Varies

For up-to-date licensing information and product-specific disclaimers, see the respective QIAGEN kit handbook or user manual. QIAGEN kit handbooks and user manuals are available at www.qiagen.com or can be requested from QIAGEN Technical Services or your local distributor.

Notes

Notes

Trademarks: QIAGEN® (QIAGEN Group).

This warranty limits our liability to replace this product in the event the product fails to perform due to any manufacturing defect. QIAGEN makes no other warranties of any kind, expressed or implied, including without limitation, warranties of merchantability or fitness for a particular purpose. QIAGEN shall not be liable for any direct, indirect, consequential or incidental damages arising out of the use, the results of use or the inability to use this product.

NOTICE TO PURCHASER

Purchase of an ELISArray Kit does not grant rights to use the kit components for reproduction of any primer pair mix, to modify kit components for resale or to use the ELISArray Kit to manufacture commercial products without written approval of QIAGEN. No other license, expressed, implied or by estoppel, is granted. U.S. patents may cover the use of certain antibodies included in the ELISArray Kit. Presently, it is not clear under U.S. laws whether commercial users must obtain licenses from the owners of the rights to these U.S. patents before using the ELISArray Kit.

Limited License Agreement

Use of these products signifies the agreement of any purchaser or user of the Single-Analyte ELISArray Kits for Rat to the following terms:

1. The Single-Analyte ELISArray Kits for Rat may be used solely in accordance with the *Single-Analyte ELISArray for Rat Handbook* and for use with components contained in the Kits only. QIAGEN grants no license under any of its intellectual property to use or incorporate the enclosed components of these Kits with any components not included within these Kits except as described in the *Single-Analyte ELISArray for Rat Handbook* and additional protocols available at www.qiagen.com.
2. Other than expressly stated licenses, QIAGEN makes no warranty that these Kits and/or their use(s) do not infringe the rights of third-parties.
3. These Kits and their components are licensed for one-time use and may not be reused, refurbished, or resold.
4. QIAGEN specifically disclaims any other licenses, expressed or implied other than those expressly stated.
5. The purchaser and user of these Kits agree not to take or permit anyone else to take any steps that could lead to or facilitate any acts prohibited above. QIAGEN may enforce the prohibitions of this Limited License Agreement in any Court, and shall recover all its investigative and Court costs, including attorney fees, in any action to enforce this Limited License Agreement or any of its intellectual property rights relating to these Kits and/or their components.

For updated license terms, see www.qiagen.com.

© 2011 QIAGEN, all rights reserved.

www.qiagen.com

Australia ■ Orders 1-800-243-800 ■ Fax 03-9840-9888 ■ Technical 1-800-243-066

Austria ■ Orders 0800-28-10-10 ■ Fax 0800-28-10-19 ■ Technical 0800-28-10-11

Belgium ■ Orders 0800-79612 ■ Fax 0800-79611 ■ Technical 0800-79556

Brazil ■ Orders 0800-557779 ■ Fax 55-11-5079-4001 ■ Technical 0800-557779

Canada ■ Orders 800-572-9613 ■ Fax 800-713-5951 ■ Technical 800-DNA-PREP (800-362-7737)

China ■ Orders 86-21-3865-3865 ■ Fax 86-21-3865-3965 ■ Technical 800-988-0325

Denmark ■ Orders 80-885945 ■ Fax 80-885944 ■ Technical 80-885942

Finland ■ Orders 0800-914416 ■ Fax 0800-914415 ■ Technical 0800-914413

France ■ Orders 01-60-920-926 ■ Fax 01-60-920-925 ■ Technical 01-60-920-930 ■ Offers 01-60-920-928

Germany ■ Orders 02103-29-12000 ■ Fax 02103-29-22000 ■ Technical 02103-29-12400

Hong Kong ■ Orders 800 933 965 ■ Fax 800 930 439 ■ Technical 800 930 425

Ireland ■ Orders 1800 555 049 ■ Fax 1800 555 048 ■ Technical 1800 555 061

Italy ■ Orders 800-789-544 ■ Fax 02-334304-826 ■ Technical 800-787980

Japan ■ Telephone 03-6890-7300 ■ Fax 03-5547-0818 ■ Technical 03-6890-7300

Korea (South) ■ Orders 080-000-7146 ■ Fax 02-2626-5703 ■ Technical 080-000-7145

Luxembourg ■ Orders 8002-2076 ■ Fax 8002-2073 ■ Technical 8002-2067

Mexico ■ Orders 01-800-7742-639 ■ Fax 01-800-1122-330 ■ Technical 01-800-7742-436

The Netherlands ■ Orders 0800-0229592 ■ Fax 0800-0229593 ■ Technical 0800-0229602

Norway ■ Orders 800-18859 ■ Fax 800-18817 ■ Technical 800-18712

Singapore ■ Orders 1800-742-4362 ■ Fax 65-6854-8184 ■ Technical 1800-742-4368

Spain ■ Orders 91-630-7050 ■ Fax 91-630-5145 ■ Technical 91-630-7050

Sweden ■ Orders 020-790282 ■ Fax 020-790582 ■ Technical 020-798328

Switzerland ■ Orders 055-254-22-11 ■ Fax 055-254-22-13 ■ Technical 055-254-22-12

UK ■ Orders 01293-422-911 ■ Fax 01293-422-922 ■ Technical 01293-422-999

USA ■ Orders 800-426-8157 ■ Fax 800-718-2056 ■ Technical 800-DNA-PREP (800-362-7737)

