

UCP HiFidelity PCR Kit

Superior, inhibitor-resistant, high-fidelity PCR for unquestionable results

Overview

Procedure

Superior
Inhibitor
Resistance

No Misleading
Background

Whole Blood
Amplification

Amplify
25–85% GC

70x > Fidelity
than Taq

Ordering
Information



Overview

Procedure

Superior Inhibitor Resistance

No Misleading Background


Whole Blood Amplification

Amplify 25–85% GC

70x > Fidelity than Taq

Ordering Information

Ultrafast procedure with simple color change confirming correct pipetting

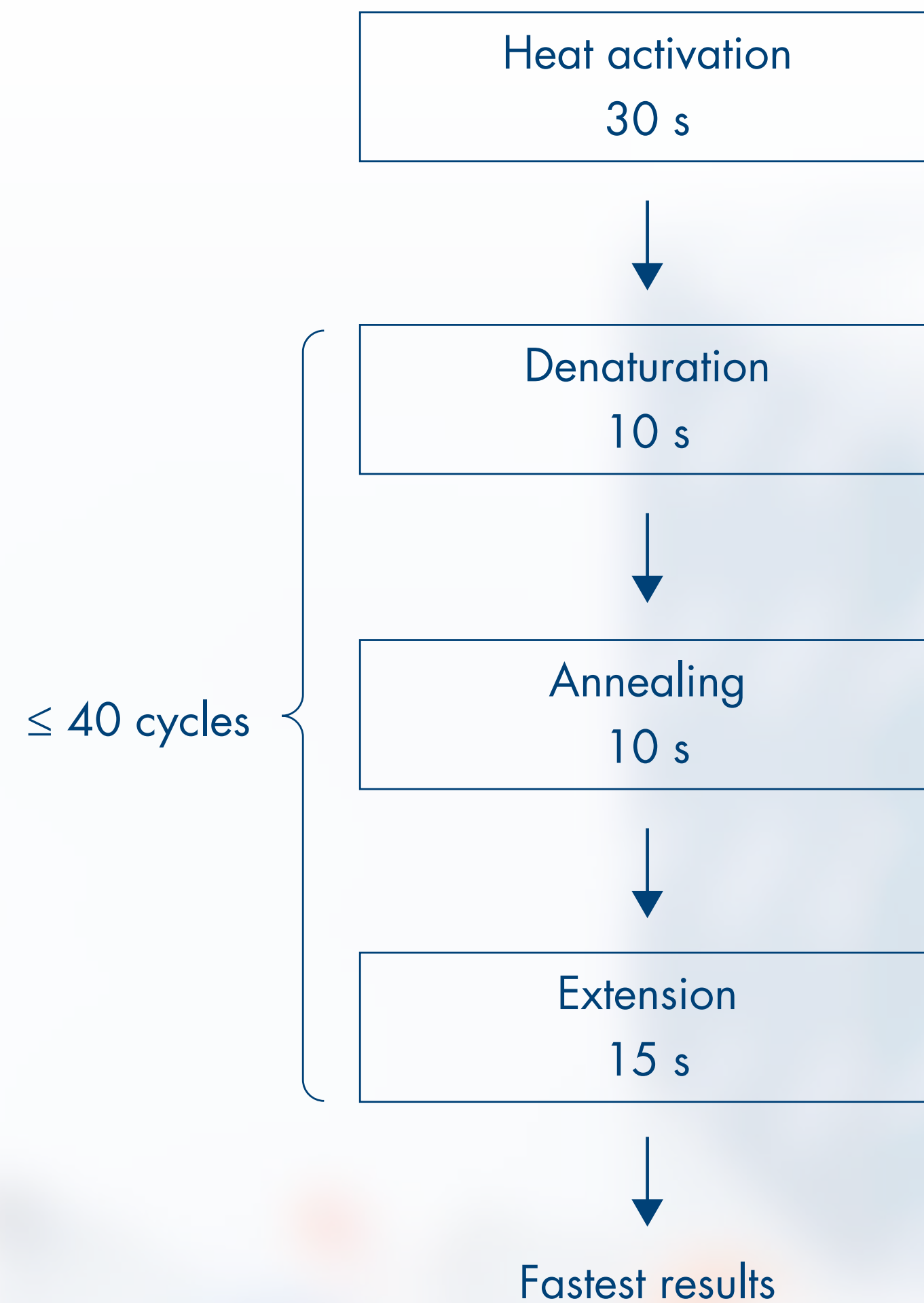

Download Quick-Start Protocol

- Unmatched convenience and speed
- Built-in pipetting control and room-temperature setup

Robust, inhibitor-resistant PCR from your most challenging samples

Overview

Procedure

Stringent Hot Start


No Misleading Background

Whole Blood Amplification

Amplify 25–85% GC

70x > Fidelity than Taq

Ordering Information

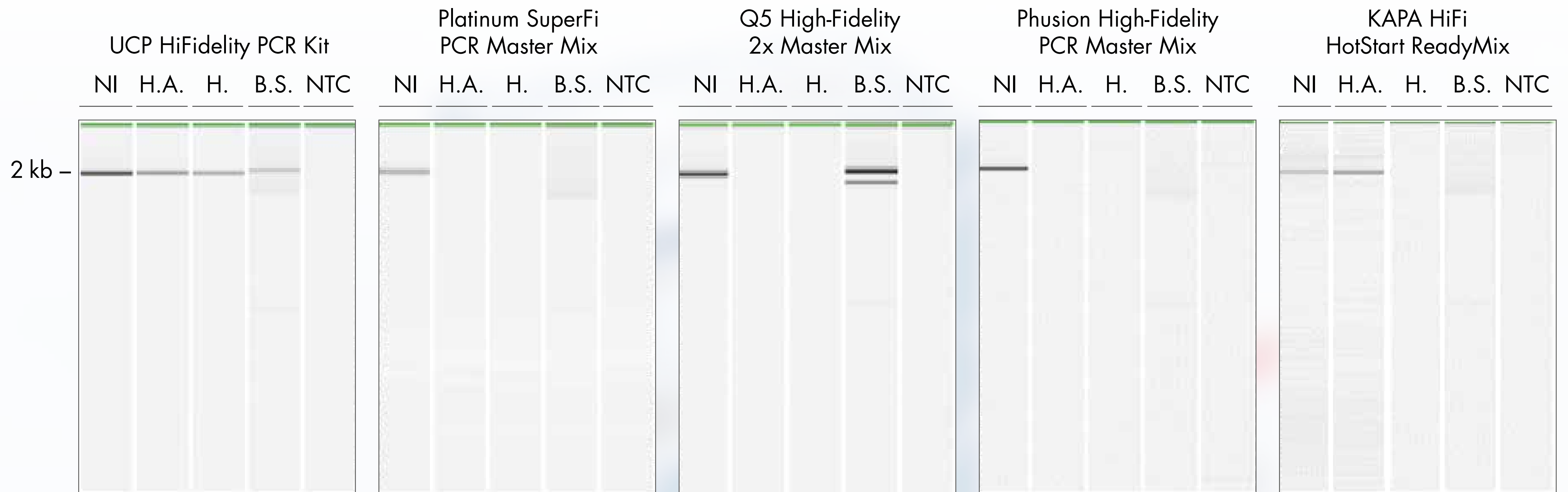


No inhibitor (N.I.), 25 ng Humic acid (H.A.), 25 μ M Heparin (H), 25 ng Bile salt (B.S.), no template control (NTC)



Superior inhibitor resistance compared to competitors

Ultra-Clean reagents for sensitive microbial detection with no background

- Overview
- Procedure
- Superior Inhibitor Resistance
- No Misleading Background
- Whole Blood Amplification
- Amplify 25–85% GC
- 70x > Fidelity than Taq
- Ordering Information

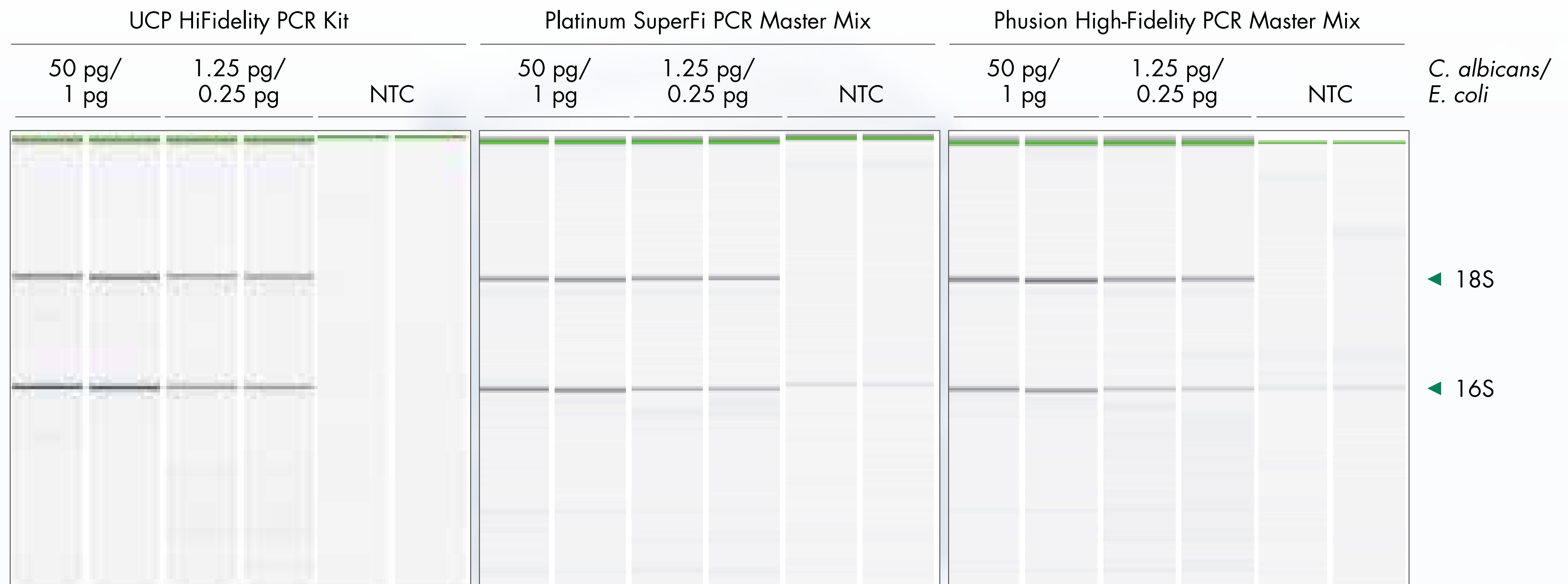

✓ No misleading background detected with Ultra-Clean reagents

Reliable amplification directly from whole blood

Overview

Procedure

Superior
Inhibitor
Resistance


No Misleading
Background

Whole Blood
Amplification

Amplify
25–85% GC

70x > Fidelity
than Taq

Ordering
Information

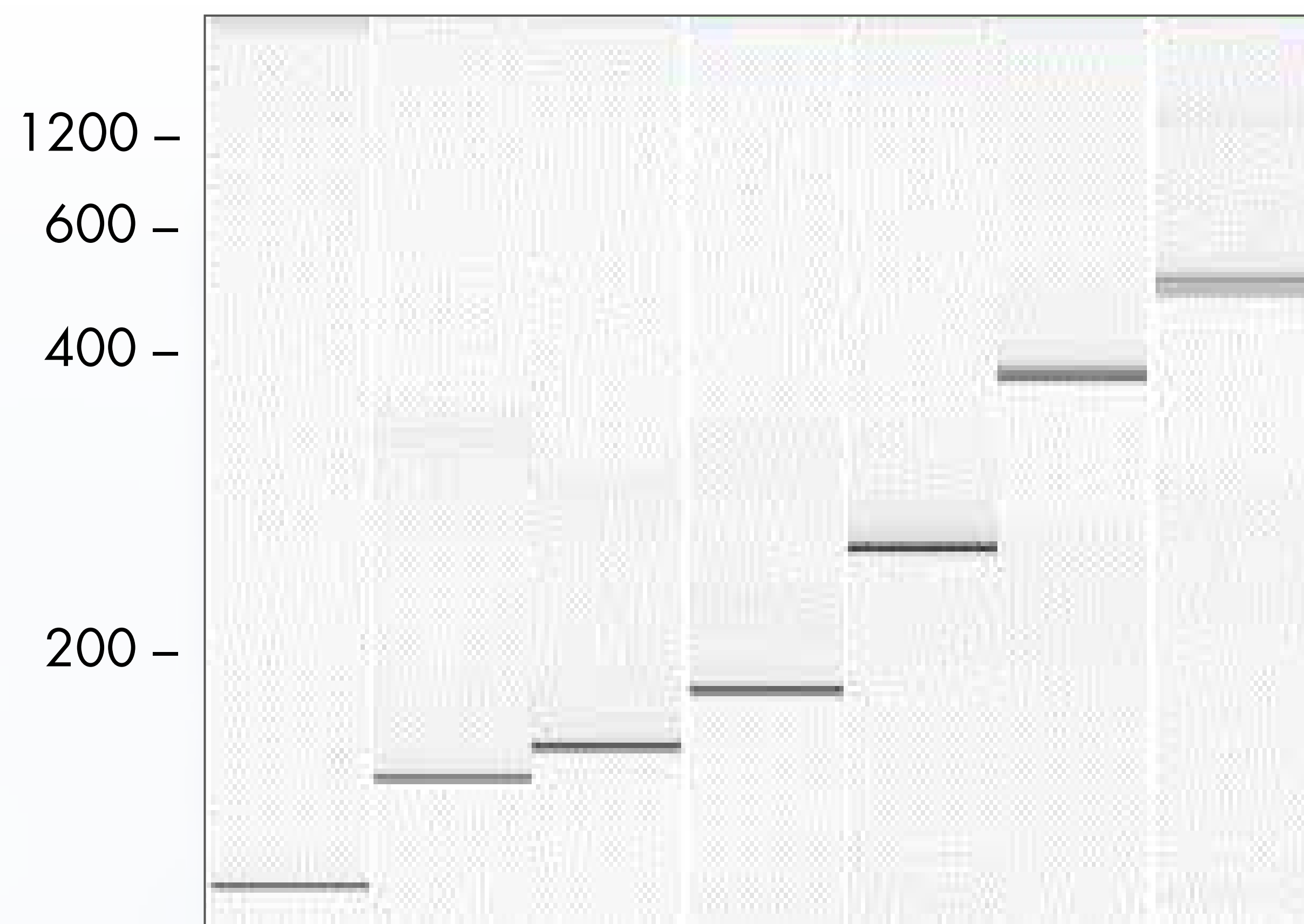

Exceptional inhibitor robustness enabling successful amplification of various targets directly from whole blood

Amplify your most challenging targets with no optimization required

Overview

Procedure

Superior Inhibitor Resistance


No Misleading Background

Whole Blood Amplification

Amplify 25–85% GC

70x > Fidelity than Taq

Ordering Information

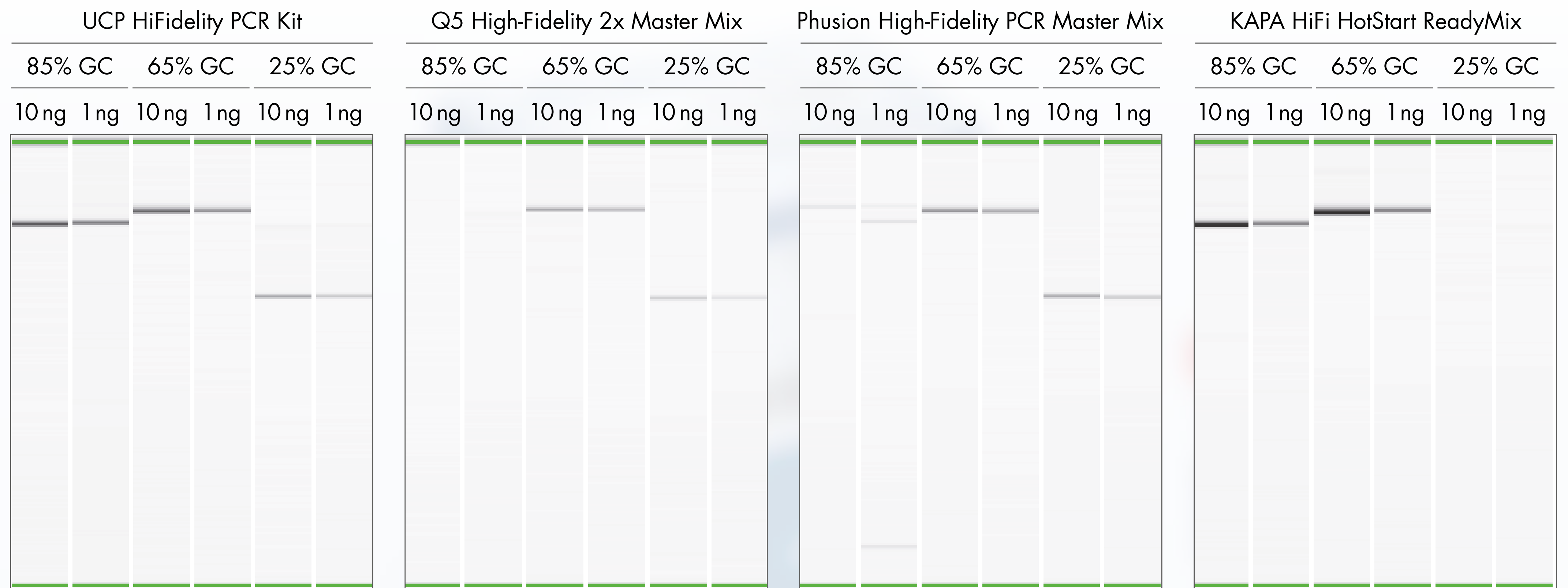

Amplify high and low GC-content templates with no protocol or buffer optimization


[Overview](#)[Procedure](#)[Superior Inhibitor Resistance](#)[No Misleading Background](#)[Whole Blood Amplification](#)[Amplify 25–85% GC](#)

70x > Fidelity than Taq

[Ordering Information](#)

Highest possible fidelity for your most rigorous applications

Fold fidelity vs. Taq

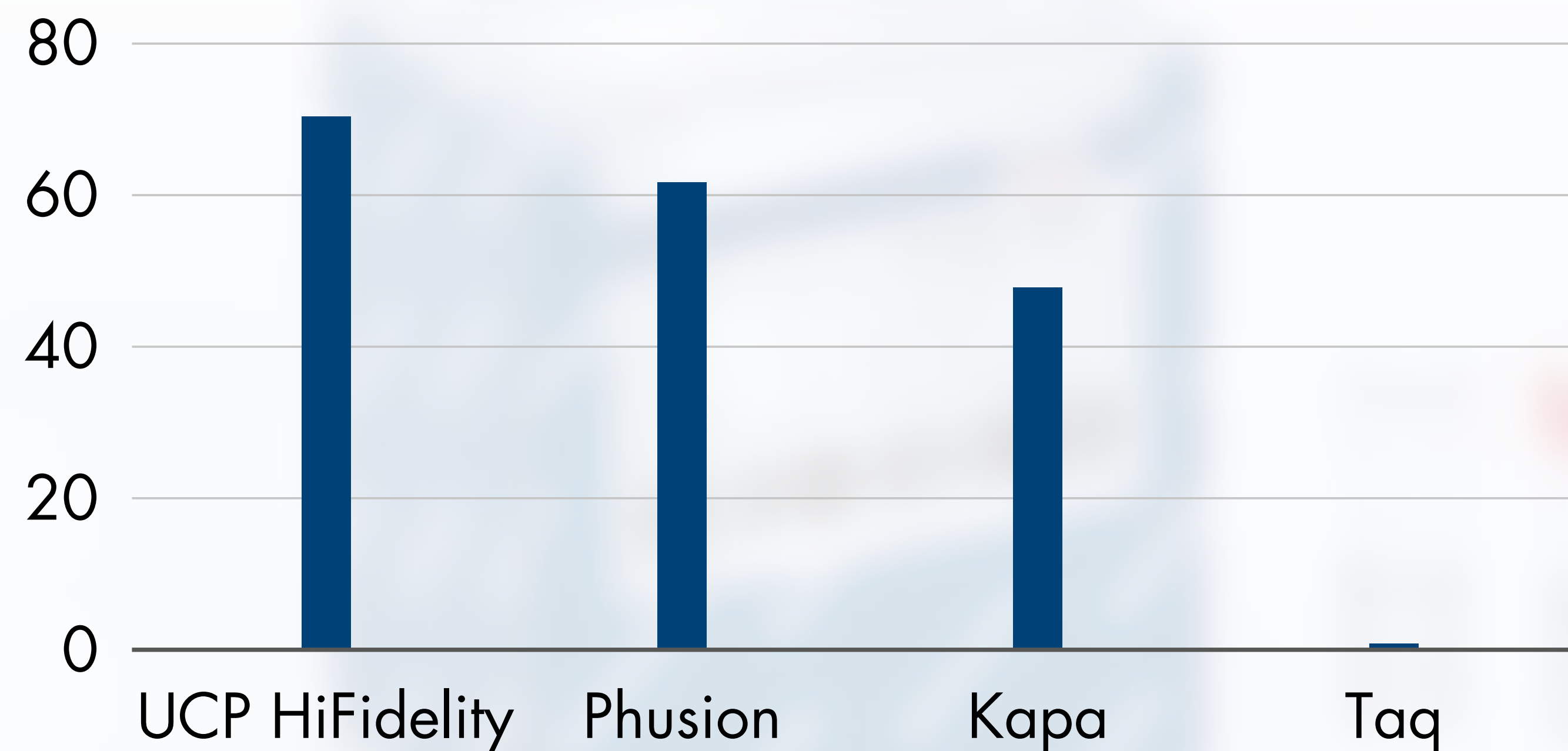

70x greater fidelity than Taq in the rigorous Lacl Fidelity test

Ordering Information

Overview
Procedure
Superior Inhibitor Resistance
No Misleading Background
Whole Blood Amplification
Amplify 25–85% GC
70x > Fidelity than Taq

Product	Contents	Cat. no.
UCP HiFidelity PCR Kit (100)	1 x 1.3 ml UCP HiFidelity MasterMix (2x), 1 x 200 µl UCP Template Tracer (25x), 1 x 50 µl UCP Master Mix Tracer (125x), 1 x 1.9 ml UCP PCR Water	202742
UCP HiFidelity PCR Kit (500)	5 x 1.3 ml UCP HiFidelity MasterMix (2x), 2 x 200 µl UCP Template Tracer (25x), 1 x 50 µl UCP Master Mix Tracer (125x), 4 x 1.9 ml UCP PCR Water	202744

For up-to-date licensing information and product-specific disclaimers, see the respective QIAGEN kit handbook or user manual. QIAGEN kit handbooks and user manuals are available at www.qiagen.com or can be requested from QIAGEN Technical Services or your local distributor.

Trademarks: QIAGEN®, Sample to Insight®, UCP™ (QIAGEN Group); Platinum®, SuperFi™, Phusion® (Thermo Fisher Scientific Inc.); Q5® (New England Biolabs).

This product is licensed from Bio-Rad Laboratories, Inc., under U.S. Pat. Nos. 6,627,424, 7,541,170, 7,560,260 and corresponding patents in other countries for use solely in DNA sequencing, DNA micro-array, and conventional PCR applications, including pre-amplification steps that are required for such applications, in the life science research and in-vitro diagnostics fields but not real-time PCR or digital PCR.

© 2018 QIAGEN, all rights reserved.

Ordering www.qiagen.com/contact | Technical Support support.qiagen.com | Website www.qiagen.com

[Visit Webpage](#)