



September 2024

## Urine Collection Procedure with PAXgene® Urine Collection Cup

Part of the PAXgene Urine Liquid Biopsy Set (cat. no. 769143)

### For collection of up to 120 mL of urine.

A minimum of 15 mL is recommended to be present in the cup to fill one tube.

The PAXgene Urine Liquid Biopsy Set is for Research Use Only. Not for use in diagnostic procedures. This product is not intended for the diagnosis, prevention, or treatment of disease.

### Preparation for specimen collection



**Step 1:** Verify you have a PAXgene Urine Collection Cup and instructions on how to collect urine.

**Step 2:** Place the collection cup aside. Wash hands thoroughly with soap and water before starting collection.

**Step 3:** Remove the blue screw cap of the PAXgene Urine Collection Cup. Place the cap on a flat surface with the "straw" facing upwards. Be careful not to touch the inside of the cap or the straw. Do not remove the warning label on top of the blue cap.

**CAUTION: SHARP NEEDLE IS LOCATED UNDER THE CAP WARNING LABEL. DO NOT REMOVE THE WARNING LABEL ON THE TOP OF THE CAP. IF THE LABEL IS REMOVED, CALL LABORATORY PROFESSIONAL TO RECEIVE A NEW CUP.**



For further details, please visit [www.preanalytix.com](http://www.preanalytix.com)


PreAnalytiX GmbH, 8634 Hombrechtikon, CH.  
The PAXgene Urine Liquid Biopsy Set and Tube are made in the UK by BD for PreAnalytiX GmbH

Unless otherwise noted, PAXgene, PreAnalytiX, the PreAnalytiX Logo and all other trademarks are property of PreAnalytiX GmbH, Hombrechtikon, CH.  
BD is trademark of Becton, Dickinson and Company  
Patent [www.preanalytix.com/patents](http://www.preanalytix.com/patents)




### Male Instructions

Prior to collection, follow instructions of the laboratory professional (for example, regarding collection time, fasting, drinking, cleaning).

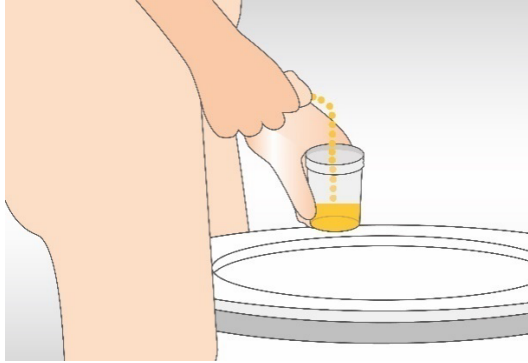

**Step 4:** Bring the collection cup closer to the end of the penis and void urine into the cup. DO NOT touch the inside or lip of the cup with the hands or any other part of the body. Once the cup is filled, void remainder of urine into the toilet.

**Important:** Please void urine only after opening the blue cap of the collection cup. Do not void into the cup through the needle hole.

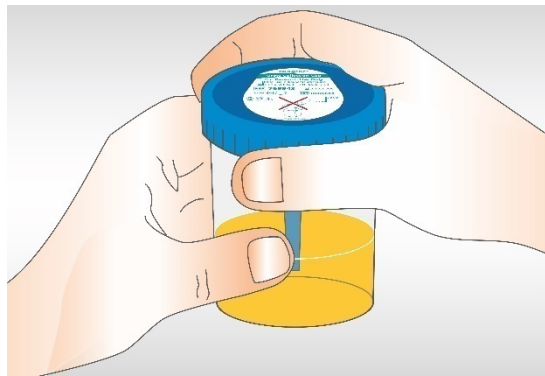

**Step 5:** Close the cup firmly with the blue cap, touching ONLY the outside surfaces of the cap and cup.


### Female Instructions

Prior to collection, follow instructions of the laboratory professional for example regarding collection time, fasting, drinking, cleaning.



**Step 4:** Bring the collection cup closer to the urinary opening and void urine into the cup. DO NOT touch the inside or lip of the cup with the hands or any other part of the body. Once the cup is filled, void remainder of urine into the toilet.

**Important:** Please void urine only after opening the blue cap of the collection cup. Do not void into the cup through the needle hole.

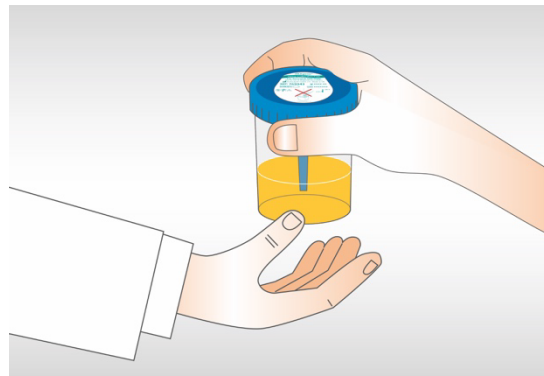

**Step 6:** Follow the instructions of the laboratory professional for immediate return of the filled urine collection cup. Keep the cup upright and avoid shaking.

For steps following collection, please refer to the *PAXgene Urine Liquid Biopsy Set Handbook* available on product webpage under **Resources** on [www.preanalytix.com](http://www.preanalytix.com) or [www.qiagen.com](http://www.qiagen.com)



For up-to-date licensing information and product-specific disclaimers, see the respective PreAnalytiX® or QIAGEN® kit handbook or user manual. PreAnalytiX and QIAGEN kit handbooks and user manuals are available at [www.preanalytix.com](http://www.preanalytix.com) and [www.qiagen.com](http://www.qiagen.com) or can be requested from QIAGEN Technical Service or your local distributor.