

Important Note

Compatible labware on the EZ2[®] Connect Fx

Dear valued customer,

We are writing to clarify what labware is compatible on the EZ2 Connect Fx with the EZ1&2[™] DNA Investigator Kit protocols. We understand the need for flexibility in your workflow and want to provide you with this information to prevent any technical issues and ensure continued success using the EZ2 Connect Fx. The following labware is compatible on the EZ2 Connect Fx using the EZ1&2 DNA Investigator Kit protocols:

Standard tip rack, sample tubes

- QIAGEN[®] Sample Tubes CB (2 mL) (QIAcube[®] accessories, cat. no. 990382)
- QIAGEN Sample Tube (2 mL) (included in EZ1&2 DNA Investigator[®] Kit, cat. no. 952034)

Standard tip rack, elution tubes

- QIAGEN Elution Tube 1.5 mL (included in EZ1&2 DNA Investigator Kit)
- Sarstedt[®] 1.5 mL elution tube (Sarstedt, cat. no. 72.692.405)
- Sarstedt 1.5 mL (Sarstedt, cat. no. 72.690.001)

Flip-cap rack, sample tubes

- QIAGEN Sample Tubes CB (2 mL) (QIAcube accessories, cat. no. 990382)
Note: May not pass optional load check due to missing lid
- QIAGEN Sample Tube (2 mL) (included in EZ1&2 DNA Investigator Kit, cat. no. 952034)
Note: May not pass optional load check due to missing lid
- Eppendorf® Safe Lock 2 mL (Eppendorf, cat. no. 0030 120.094)
- Sarstedt SafeSeal 2 mL (Sarstedt, cat. no. 72.695.500)
- Sarstedt 2 mL sample tube (Sarstedt, cat. no. 72-694.406)
Note: May not pass optional load check due to missing lid
- QIAGEN Sample Tube 2 mL (included in Investigator Lyse&Spin Basket Kit, cat nos. 19597 and 19598)

Note: Ensure that flip-cap tubes are sitting flush after inserting the lids.

Note: We recommend not using the 1.5 mL QIAcube Connect elution tube in the sample tube position of the EZ2 Connect Fx, unless using protocols dedicated to these tubes.

Flip-cap rack, elution tubes

- QIAGEN Elution Tube 1.5 mL (included in EZ1&2 DNA Investigator Kit, cat. no. 953024)
Note: May not pass optional load check due to missing lid
- Sarstedt 1.5 mL EZ1® screw-cap elution tube (Sarstedt, cat. no. 72.692.405)
Note: May not pass optional load check due to missing lid
- Sarstedt 1.5 mL (Sarstedt, cat. no. 72.690.001)

- Eppendorf 1.5 mL 3810X (Eppendorf, cat. no. 0030 125.150)
- Eppendorf Safe Lock 1.5 mL (Eppendorf, cat. no. 0030 120.086)

Large volume rack, sample tubes

- Sarstedt screw cap tube, working volume: 7 mL, 82 x 13 mm (Sarstedt, cat. no. 60.550.109)
- QIAGEN 7 mL Large-Volume Tubes (cat. no. 951954)

Large volume rack, elution tubes

- QIAGEN Elution Tube 1.5 mL (included in EZ1&2 DNA Investigator Kit, cat. no. 952034)
- Sarstedt 1.5 mL EZ1 screw-cap elution tube (Sarstedt, cat. no. 72.692.405)
- Sarstedt 1.5 mL (Sarstedt, cat. no. 72.690.001)

Heater position

- Sarstedt 2 mL tube (cat. no. 72.693.005)
- QIAGEN Sample Tubes CB (2 ml) (QIAcube accessories, cat. no. 990382)

Note: QIAGEN Sample Tube 2.0 ml (skirted) that is included in EZ1&2 DNA Investigator Kit is not compatible in heater position of the EZ2 Connect Fx.

If you have further questions, please contact your local QIAGEN representative or contact our Technical Support Center at www.qiagen.com

Best regards,

QIAGEN

Trademarks: QIAGEN®, Sample to Insight®, QIAcube®, QIAxcel®, EZ1®, EZ2®, EZ1&2™, Investigator® (QIAGEN Group); Eppendorf® (Eppendorf AG); Sarstedt® (Sarstedt AG and Co.). Registered names, trademarks, etc. used in this document, even when not specifically marked as such, are not to be considered unprotected by law.

01/2024 HB-3546-001 © 2024 QIAGEN, all rights reserved.