TissueLyser III Safety Instructions and Quick-Start Guide

Contents

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EN

Introduction

Thank you for choosing the TissueLyser III system.

Before using the TissueLyser III, it is essential you read this quick-start guide carefully and pay particular attention to the safety information. The instructions and safety information in this guide must be followed to ensure safe operation of the TissueLyser III and to maintain the TissueLyser III in a safe condition. For more detailed information and instructions, visit the product page of the TissueLyser III on [qiagen.com](https://www.qiagen.com) and download the *TissueLyser III User Manual*.

Safety Information

This instrument is intended for professional use.

The TissueLyser III is to be used only by personnel specifically instructed and trained in biology laboratory techniques and familiar with this technology.

Before using the TissueLyser III, it is essential that you read this quick-start guide carefully. The instructions and safety information in the quick-start guide must be followed to ensure safe operation of the instrument and to maintain the instrument in a safe condition.

The following types of safety information appear throughout the *TissueLyser III Safety Instructions and Quick Start Guide*.

<p>WARNING</p> 	<p>The term WARNING is used to inform you about situations that could result in personal injury to you or others.</p> <p>Details about these circumstances are given in a box like this one.</p>
<p>CAUTION</p> 	<p>The term CAUTION is used to inform you about situations that could result in damage to an instrument or other equipment.</p> <p>Details about these circumstances are given in a box like this one.</p>

The guidance provided in this guide is intended to supplement, not supersede, the normal safety requirements prevailing in the user's country.

Please be aware that you may be required to consult your local regulations for reporting serious incidents that have occurred in relation to the device to the manufacturer and/or its authorized representative (only applicable for CE-marked devices with an authorized representative established in the EU) and the regulatory authority in which the user and/or the patient is established.

Proper use

<p>WARNING</p> 	<p>Risk of personal injury and material damage</p> <p>Improper use of the TissueLyser III may cause personal injuries or damage to the instrument. Do not attempt to move the TissueLyser III during operation.</p> <p>The TissueLyser III should only be operated by qualified personnel who have been appropriately trained.</p> <p>Servicing of the TissueLyser III should only be performed by QIAGEN® or an authorized agent (service technicians). Do not carry out any modifications of the instrument.</p> <p>QIAGEN shall not be held responsible for any injuries or damages resulting from nonobservance of the instructions and safety information in this guide.</p>
<p>WARNING</p>  	<p>The sound level may be high depending on the type of material, the number of balls used, the set grinding frequency, and the grinding time. Excess noise in terms of intensity and duration can lead to impairments or permanent damage to hearing. Wear hearing protection.</p> <p>Excess noise can also lead to acoustic warning signals of other instruments not being heard. Add visual signals if necessary.</p>
<p>CAUTION</p> 	<p>Damage to the instrument</p> <p>Use of adapters and plasticware other than those described in the <i>TissueLyser III User Manual</i> may lead to damage to the clamps of the TissueLyser III.</p>

CAUTION**Damage to the instrument**

Avoid spilling water or chemicals onto the Tissuelyser III. Damage caused by water or chemical spillage will void your warranty.

CAUTION**Damage to the instrument**

Operation of the Tissuelyser III with the transport security device installed, or transport of the Tissuelyser III without the transport security device installed, can lead to damage to the mechanical components of the instrument.

CAUTION**Damage to the instrument**

Operation of the Tissuelyser III at a temperature and humidity outside the range specified may lead to damage to its electronic and mechanical components as well as a reduction in instrument performance.

WARNING**Risk of personal injury and material damage through falling of the instrument**

Improper handling of the Tissuelyser III during transport and installation can lead to user injury. Particularly when lifting the instrument above head height, the instrument can fall and lead to serious injury.

Do not transport the instrument alone. Transport the instrument close to the floor and avoid lifting above the head.

Install the instrument on a sufficiently large, solid, and stable surface that is protected against oscillations. Ensure that all instrument feet are sitting stably and evenly on the surface.

Electrical safety

<p>WARNING</p> 	<p>Electrical hazard</p> <p>Any interruption of the protective conductor (earth/ground lead) inside or outside the Tissuelyser III or disconnection of the protective conductor terminal is likely to make the Tissuelyser III dangerous.</p> <p>Intentional interruption is prohibited.</p> <p>Take note of the values for voltage and frequency on the Tissuelyser III type plate and only connect the Tissuelyser III to a fitting power grid.</p> <p>Only connect the Tissuelyser III to the power grid with the enclosed power cable and to a power outlet with protective conductor (earth/ground).</p> <p>Before each usage, check the power cord and the plug for damage. Never use the instrument with damaged power cord or plug.</p> <p>Lethal voltages inside the Tissuelyser III.</p> <p>When the Tissuelyser III is connected to line power, terminals may be live and opening the Tissuelyser III or removing parts is likely to expose live parts.</p> <p>Install the instrument in a way that allows easy access to the power cord connector and the power switch.</p>
---	--

To ensure satisfactory and safe operation of the Tissuelyser III, follow the advice below:

- The line power cord must be connected to a line power outlet that has a protective conductor (earth/ground).
- Do not adjust or replace internal parts of the Tissuelyser III.
- Do not operate the Tissuelyser III with any covers or parts removed.
- If liquid has spilled inside the Tissuelyser III, switch off the Tissuelyser III, disconnect it from the power outlet, and contact QIAGEN Technical Services.

- If the TissueLyser III becomes electrically unsafe, prevent other personnel from operating it, and contact QIAGEN Technical Services; the TissueLyser III may be electrically unsafe when:
 - It or the line power cord appears to be damaged.
 - It has been stored under unfavorable conditions for a prolonged period.
 - It has been subjected to severe transport stresses.

If the TissueLyser III is not in use, switch it off at the power switch. If the TissueLyser III will not be used for a prolonged period of time, unplug the power cord from the power outlet.

Environment

<p>WARNING</p> 	<p>Explosive atmosphere</p> <p>The TissueLyser III is not designed for use in an explosive atmosphere.</p>
<p>WARNING</p> 	<p>Risk of explosion</p> <p>The TissueLyser III is intended for use with reagents and substances supplied with QIAGEN kits. Use of other reagents and substances may lead to fire or explosion.</p>
<p>CAUTION</p> 	<p>Damage to the instrument</p> <p>Direct sunlight may bleach parts of the instrument and cause damage to plastic parts. The TissueLyser III must be located out of direct sunlight.</p>

Biological safety

Specimens and reagents containing materials from humans or animals should be treated as potentially infectious. Use safe laboratory procedures as outlined in publications such as *Biosafety in Microbiological and Biomedical Laboratories*, CDC (1).

Samples

Samples may contain infectious agents. You should be aware of the health hazard presented by such agents and should use, store, and dispose of such samples according to the required safety regulations.

<p>WARNING</p> 	<p>Samples containing infectious agents</p> <p>Some samples used with the TissueLyser III may contain infectious agents. Handle such samples with the greatest of care and in accordance with the required safety regulations.</p> <p>Always wear safety glasses, 2 pairs of gloves, and a lab coat.</p> <p>The responsible body (e.g., laboratory manager) must take the necessary precautions to ensure that the surrounding workplace is safe and that TissueLyser III operators are suitably trained and not exposed to hazardous levels of infectious agents as defined in the applicable Safety Data Sheets (SDSs) or OSHA,* ACGIH,† or COSHH‡ documents.</p> <p>Venting for fumes and disposal of wastes must be in accordance with all national, state, and local health and safety regulations and laws.</p>
---	--

* OSHA: Occupational Safety and Health Administration (United States of America).

† ACGIH: American Conference of Government Industrial Hygienists (United States of America).

‡ COSHH: Control of Substances Hazardous to Health (United Kingdom).

Chemicals

<p>WARNING</p> 	<p>Hazardous chemicals</p> <p>Some chemicals used with the TissueLyser III may be hazardous or may become hazardous after completion of the sample disruption process.</p> <p>Always wear safety glasses, gloves, and a lab coat.</p> <p>The responsible body (e.g., laboratory manager) must take the necessary precautions to ensure that the surrounding workplace is safe and that TissueLyser III operators are not exposed to hazardous levels of toxic substances (chemical or biological) as defined in the applicable Safety Data Sheets (SDSs) or OSHA,[*] ACGIH,[†] or COSHH[‡] documents.</p> <p>Venting for fumes and disposal of wastes must be in accordance with all national, state, and local health and safety regulations and laws.</p>
---	---

* OSHA: Occupational Safety and Health Administration (United States of America).

† ACGIH: American Conference of Government Industrial Hygienists (United States of America).

‡ COSHH: Control of Substances Hazardous to Health (United Kingdom).

WARNING**Liquid nitrogen**

Works with liquid nitrogen pose risks due to oxygen displacement and cryogenic burns. Ensure proper ventilation of the room and air monitoring when handling or storing liquid nitrogen. Avoid spills by always keeping vessels upright and do not touch with bare skin vessels that are or have been in exposed to liquid nitrogen. Wear appropriate protective equipment (safety glasses, gloves, lab coat, and safety shoes).

The low temperature of liquid nitrogen can change the mechanical properties of materials. Nitrogen-cooled plastic vessels can become brittle and break when used in the TissueLyser III. Only use liquid nitrogen with vessels that have been cleared for such use in the TissueLyser III.

The responsible body (e.g., laboratory manager) must take the necessary precautions to ensure that the surrounding workplace is safe and that TissueLyser III operators are not exposed to the dangers posed by liquid nitrogen as defined in the applicable Safety Data Sheets (SDSs) or OSHA,* ACGIH,[†] or COSHH[‡] documents.

* OSHA: Occupational Safety and Health Administration (United States of America).

[†] ACGIH: American Conference of Government Industrial Hygienists (United States of America).

[‡] COSHH: Control of Substances Hazardous to Health (United Kingdom).

Toxic fumes

If working with volatile solvents, toxic substances, etc., you must provide an efficient laboratory ventilation system to remove vapors that may be produced.

WARNING**Toxic fumes**

Do not use bleach to clean or disinfect used labware. Bleach in contact with salts from the buffers used can produce toxic fumes.

Waste disposal

Used labware, such as sample tubes, may contain hazardous chemicals or infectious agents from the sample disruption process. Such wastes must be collected and disposed of properly according to local safety regulations. For information on how to dispose of the Tissuelyser III, see Appendix A of the *Tissuelyser III User Manual*.

Mechanical and thermal hazards

WARNING 	Moving parts Do not leave the Tissuelyser III unattended during operation.
---	--

To ensure safety, the Tissuelyser III will not operate until the hood is closed.

WARNING 	Moving parts Damage or personal injury can occur if the Tissuelyser Adapter Set or Grinding Jar Set is not securely attached. Ensure that both adapters and grinding jars are of similar weight. Do not operate the Tissuelyser III with only a single adapter or grinding jar installed.
---	--

WARNING 	Danger of burns and scalds Depending on the sample and the disruption parameters, grinding jars and samples can heat up considerably during the disruption process. Only touch the grinding jars with protective gloves and wait for equilibration to room temperature before opening the jars.
---	---

Maintenance safety

<p>WARNING</p> 	<p>Risk of electric shock</p> <p>Do not open any panels on the TissueLyser III. Only perform cleaning with liquids after disconnecting the instrument from the power grid. Use liquids only on cloth, never free-flowing.</p> <p>Risk of personal injury and material damage</p> <p>Only perform maintenance that is specifically described in the <i>TissueLyser III User Manual</i>.</p>
<p>CAUTION</p> 	<p>Damage to the instrument</p> <p>Do not use bleach, solvents, or reagents containing acids, alkalis, or abrasives to clean the TissueLyser III.</p> <p>Avoid spilling water or chemicals onto the TissueLyser III. Damage caused by water or chemical spillage will void your warranty.</p>

Symbols on the TissueLyser III

Symbol	Location	Description
	<p>Type plate on the back of the instrument</p>	<p>WEEE about the disposal of waste electrical and electronic equipment for Europe and rest of the world.</p>
	<p>Type plate on the back of the instrument</p>	<p>Legal manufacturer</p>
	<p>Type plate on the back of the instrument</p>	<p>Date of manufacture</p>

Symbol	Location	Description
	Type plate on the back of the instrument	Consult instructions for use.
	Type plate on the back of the instrument	CE mark for Europe
	Type plate on the back of the instrument	UKCA mark for the United Kingdom
	Type plate on the back of the instrument	Mark of the TÜV Rheinland
	Type plate on the back of the instrument	RCM (former C-Tick) for Australia (supplier identification N17965)
	Type plate on the back of the instrument	RoHS mark for China (the restriction of the use of certain hazardous substances in electrical and electronic equipment)
	Type plate on the back of the instrument	Instrument serial number
	Type plate on the back of the instrument	Global trade item number
	Type plate on the back of the instrument	Unique device identifier

Symbol	Location	Description
	Type plate on the back of the instrument	QIAGEN catalogue number
	Type plate on the back of the instrument	Label number
	Type plate on the back of the instrument	See warnings and precautions
	On the back of the instrument	Wear hearing protection
	On the back of the instrument	Refer to user manual

Installation Procedures

Unpacking the Tissuelyser III

Before unpacking the Tissuelyser III, check whether the cardboard container is damaged. In case of damage, contact the transporter of the container. The Tissuelyser III must not be knocked, shaken, or thrown during transport; otherwise, its electronic and mechanical components may be damaged.

Make a note of how the Tissuelyser III is packed. If returning the instrument to QIAGEN for repair, it needs to be repacked using the original packaging. Lift the Tissuelyser III with two people, placing both hands below the base of the instrument.

After unpacking the TissueLyser III, check that the following items are supplied:

- TissueLyser III
- Power cords for different countries
- *TissueLyser III Safety Instructions and Quick-Start Guide*
- Tool to remove the transport security device

If anything is missing, contact QIAGEN Technical Services.

Check that the TissueLyser III is not damaged. If anything is damaged, contact QIAGEN Technical Services. Make sure that the TissueLyser III has equilibrated to ambient temperature before operating it. The TissueLyser III should be stored dry when not in use.

Retain the original packaging in case you need to return the TissueLyser III to QIAGEN for repair. Using the original packaging minimizes damage during transportation of the TissueLyser III and avoids invalidation of your warranty.

Installation requirements

Removal of the transport security devices

To protect the TissueLyser III during transport, transport security devices are located on the underside of the instrument and around the clamps. It is important to remove the transport security devices before usage of the TissueLyser III to avoid damage to the instrument. The required tool is delivered together with the TissueLyser III.

To remove the transport security devices, follow the following steps:

- On the underside of the instrument, an arrow marks the location of the transport security device (Figure 1).

- Rotate the Tissuelyser III to the side and place it on a soft surface. Ensure that the instrument is protected against tilting at all times.
- Use the tool to remove the transport security device. It is a screw located in a recess at the center of the baseplate.
- Rotate the Tissuelyser III back onto its feet and remove the foam block fixing the clamps (Figure 2).
- Carefully store the transport security devices for future transport of the Tissuelyser III.

Location of the transport security device

Figure 1. Bottom transport security device. The first transport security device at the underside of the Tissuelyser III is comprised of a screw that has to be removed with the tool delivered with the instrument. Ensure that the transport security device is removed as depicted in the figure before operating the Tissuelyser III.

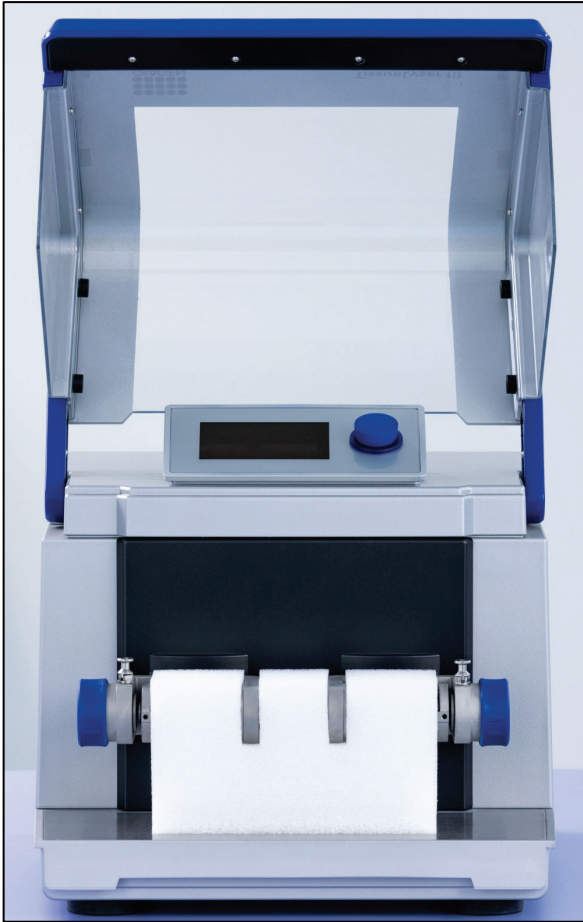
Figure 2 Inner transport security device. The second transport security device comprises of a foam block between the clamps. Remove the block.

Site requirements

Place the Tissuelyser III on a flat, stable surface, and ensure that there is sufficient space around and under the instrument for ventilation. The Tissuelyser III is for indoor use only under the following conditions:

- Ambient temperature within the range of 5–40°C (41–104°F)
- Up to 2000 m (6500 ft.) above mean sea level
- Maximum relative humidity of 80% for temperatures up to 31°C, linearly decreasing down to a relative humidity of 50% at 40°C

Power requirements

The power line to the Tissuelyser III should be voltage regulated and surge protected. Make sure that the voltage rating of the Tissuelyser III (see the plate on the back of the instrument) is compatible with the AC voltage available at the installation site.

Grounding requirements

To protect operating personnel, the National Electrical Manufacturers Association recommends that the Tissuelyser III be correctly grounded (earthed). The instrument should be plugged into an AC power outlet that has a ground (earth) connection.

Reference

1. Meechan PJ, Hatcher B, Potts J. *Biosafety in Microbiological and Biomedical Laboratories (BMBL)*. 6th ed. Centers for Disease Control and Prevention; 2020.
<https://www.cdc.gov/labs/BMBL.html>

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Document Revision History

Revision	Description
May 2023	Initial release

For up-to-date licensing information and product-specific disclaimers, see the respective QIAGEN kit handbook or user manual. QIAGEN kit handbooks and user manuals are available at www.qiagen.com or can be requested from QIAGEN Technical Services or your local distributor.

This page intentionally left blank

Trademarks: QIAGEN®, Sample to Insight® (QIAGEN Group). Registered names, trademarks, etc. used in this document, even when not specifically marked as such, are not to be considered unprotected by law.

May-2023 HB-3240-001 1 130034 © 2023 QIAGEN, all rights reserved.