

QuantiNova[®] SYBR[®] Green PCR Kit

The QuantiNova SYBR Green PCR Kit (cat. nos. 208052, 208054, 208056) should be stored immediately upon receipt at -30°C to -15°C in a constant-temperature freezer and protected from light. QuantiNova SYBR Green PCR Master Mix can also be stored protected from light at $2-8^{\circ}\text{C}$ for up to 6 months, depending on the expiration date.

Further information

- *QuantiNova SYBR Green PCR Handbook*: www.qiagen.com/HB-1698
- Safety Data Sheets: www.qiagen.com/safety
- Technical assistance: support.qiagen.com

Notes before starting

- This protocol is optimized for quantification of gDNA or cDNA targets using SYBR Green I with any real-time cyclers and condition for fluorescence normalization. ROX[™] dye is required for various cyclers at the following concentrations:
 - No requirement for ROX dye: Rotor-Gene[®], Bio-Rad[®] CFX, Roche LightCycler[®] 480 and Agilent[®] Technologies Mx instruments
 - Low concentration of ROX dye: Applied Biosystems[®] 7500, ViiA[®]7 and QuantStudio[®] Real-Time PCR Systems
 - High concentration of ROX dye: ABI PRISM[®] 7000, Applied Biosystems 7300, 7900 and StepOne[™] Real-Time PCR Systems
- ROX Reference Dye is provided as a separate tube of passive reference dye for normalization of fluorescent signals on all real-time cyclers from Applied Biosystems. It should be diluted 1:10 in the 1x reaction when used on instruments requiring a high ROX concentration and 1:200 for instruments requiring a low ROX concentration.
- The dye in QuantiNova Yellow Template Dilution Buffer allows tracking of pipetted samples in the qPCR. When the template is added to the blue QuantiNova SYBR Green PCR Master Mix, the color changes from blue to green, indicating the successful inclusion

of template. The use of this buffer is optional. It is provided as 100x concentrate and should be diluted (using water or Tris buffer) to obtain a final concentration of 1x within samples. To generate a template dilution series (e.g., for absolute quantification or determination of PCR efficiency), dilute the 100x concentrate (using water or Tris buffer) to obtain a final concentration of 1x QuantiNova Yellow Template Dilution Buffer. The buffer does not affect the sample stability and qPCR.

- For the highest efficiency in real-time PCR using SYBR Green I, amplicons should ideally be 60–200 bp in length. For amplicons >200 bp, a longer annealing/extension of 30 s can improve results.
- The two-step PCR cycling protocol, which has a denaturation step at 95°C and a combined annealing/extension step at 60°C, will also work for primers with a T_m well below 60°C.
- Always start with the cycling conditions and primer concentrations specified in this protocol.
- The PCR must start with an initial incubation step of 2 min at 95°C to activate the QuantiNova DNA Polymerase.

Procedure

1. Thaw QuantiNova SYBR Green PCR Master Mix, QuantiNova Yellow Template Dilution Buffer, template gDNA or cDNA, primers, probes, ROX Reference Dye (if required), and RNase-free water. Mix the individual solutions.
2. Prepare a reaction mix according to Table 1. Due to the hot start, it is not necessary to keep samples on ice during reaction setup or while programming the real-time cycler.
3. Mix the reaction mix thoroughly, and dispense appropriate volumes into PCR tubes or the wells of a PCR plate.

Table 1. Reaction setup

Component	Volume/reaction (µL)		
	96-well block, Rotor-Gene	384-well block	Final Concentration
Reaction mix 2x SYBR Green PCR Master Mix	10	5	1x
ROX Reference Dye (Applied Biosystems instruments only)	2/0.1*	1/0.05 *	1x
Primer A [†]	Variable	Variable	1 µL
Primer B [†]	Variable	Variable	Variable
RNase-free water	Variable	Variable	–
Template gDNA or cDNA (added at step 4)	Variable	Variable	≤100 ng/reaction
Total reaction volume	20	10	15 µL

* Corresponds to a 1:10 dilution for high ROX cyclers (i.e., ABI PRISM 7000, Applied Biosystems 7300, 7900 and StepOne Real-Time PCR Systems and a 1:200 dilution for low-ROX dye cyclers (i.e., Applied Biosystems 7500 and ViiA7 Real-Time PCR Systems).

[†] If using QuantiNova LNA Assays, the final concentration in the reaction should be 1x.

4. Add template gDNA or cDNA (≤100 ng/reaction) to the individual PCR tubes or wells containing the reaction mix.

Note: For two-step RT-PCR, the volume of the cDNA added (from the undiluted reverse transcription reaction) should not exceed 10% of the final PCR volume.

5. Program the real-time cyclers according to Table 2.

Note: Data acquisition should be performed during the combined annealing/extension step.

6. Place the PCR tubes or plates in the real-time cyclers and start the cycling program.

Table 2. Cycling conditions

Step	Time	Temperature (°C)	Ramp rate
PCR initial heat activation	2 min	95	1 μ L
2-step cycling:			
Denaturation	5 s	95	Maximal/fast mode
Combined annealing/extension	10 s*	60 [†]	Maximal/fast mode
Number of cycles	35–40 [‡]		
Melting curve analysis [§]			

* For amplicons >200 bp a longer annealing/extension of 30 s is recommended. If your cyclers does not accept this short time for data acquisition, use the shortest acceptable time.

[†] This temperature should also be used for QuantiNova LNA Assays and for all primer sets with a T_m well below 60°C.

[‡] The number of cycles depends on the amount of template DNA.

[§] Melting curve analysis is an analysis step

7. Perform a melting curve analysis of the PCR product(s).

Note: We strongly recommend routinely performing this analysis, which is built into the software of real-time cyclers, to verify the specificity and identity of PCR products.

Optional: Check the specificity of PCR product(s) by agarose gel electrophoresis.

Document Revision History

Date	Changes
04/2024	Converted to the new branding template. Replaced QuantiTect Primer Assays with QuantiNova LNA Assays.



Scan QR code for *QuantiNova SYBR Green PCR handbook*.

For up-to-date licensing information and product-specific disclaimers, see the respective QIAGEN kit handbook or user manual.

Trademarks: QIAGEN®, Sample to Insight®, QuantiNova®, QuantiTect®, Rotor-Gene® (QIAGEN Group); ROX™, ViiA7®, ABI PRISM®, SYBR®, StepOne™ (Life Technologies Corporation); Bio-Rad® (Bio-Rad Laboratories, Inc), LightCycler® (Roche Group); Agilent® (Agilent Technologies); Applied Biosystems® (Applied Biosystems Corporation or its subsidiaries). Registered names, trademarks, etc. used in this document, even when not specifically marked as such, are not to be considered unprotected by law. This product is provided under an intellectual property license from Life Technologies Corporation. The sale of this product is expressly conditioned on the buyer using the product or its components only in research conducted by the buyer (whether the buyer is an academic or for-profit entity). For information on purchasing a license to this product for purposes other than research, contact Life Technologies Corporation, 5791 Van Allen Way, Carlsbad, CA 92008 USA or Qullicensing@lifetechn.com.

1135228 04/2024 HB-1669-003 © 2024 QIAGEN, all rights reserved.