

# EVALUATION OF THE PAXGENE® TISSUE SYSTEM: PRESERVATION OF MORPHOLOGY AND GENE EXPRESSION IN HUMAN MELANOMA

Peter Hesse, Ralf Wyrich, Thorsten Voss, Lynne Rainen and Daniel Grözl,


PreAnalytiX GmbH, Hombrechtikon, CH

## Introduction


### PAXgene Tissue System

- Formalin-free tissue fixation method
- For simultaneous preservation of tissue morphology and biomolecules
- Allows analysis of biomarkers from the same sample used for histopathology
- Two reagent system for tissue fixation and stabilization
- Consists of the PAXgene Tissue Container, PAXgene Tissue Kits and supplementary protocols (www.preanalytix.com)
- Currently under evaluation by the EU FP7 funded project SPIDIA

### PAXgene Tissue Workflow

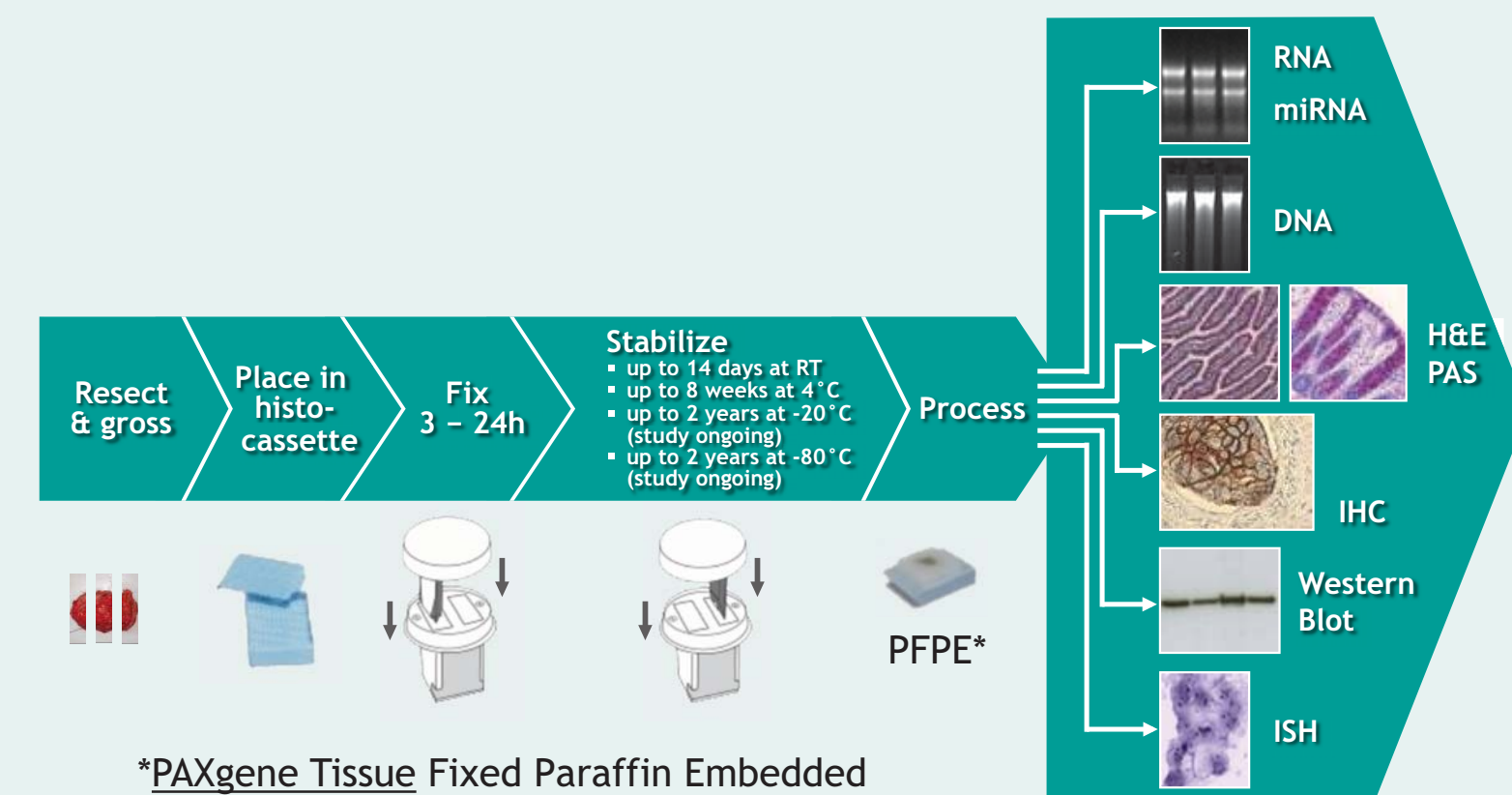


## Materials and Methods

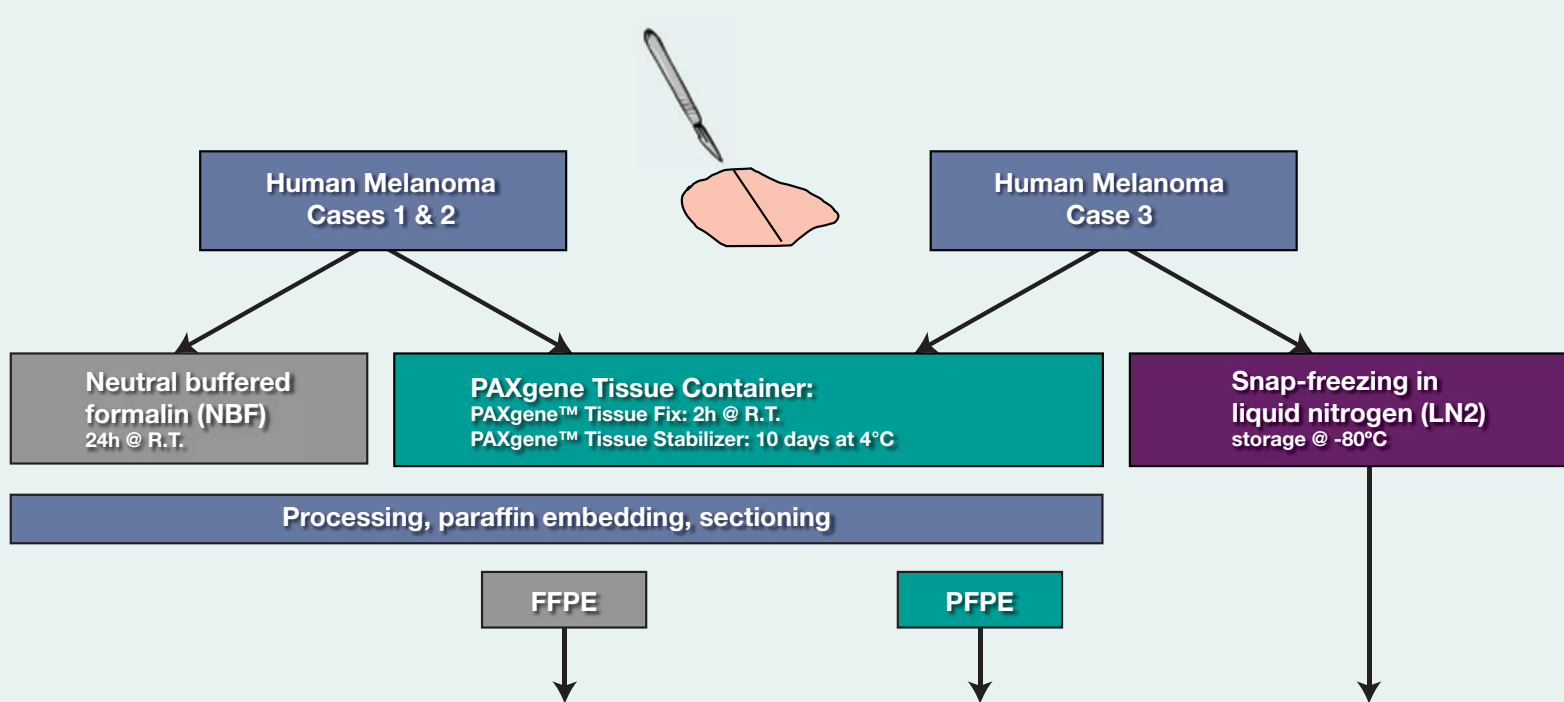

Staining	Hematoxylin & Eosin, Immunohistochemical staining for Ki-67 and S100	-
RNA purification	RNeasy® FFPE (QIAGEN)	PAXgene Tissue RNA (PreAnalytiX)
RNA incl. miRNA purification	miRNeasy® FFPE (QIAGEN)	PAXgene Tissue miRNA (PreAnalytiX)
RT-qPCR array	TaqMan® Array Gene Signature 96-Well Plate human endogenous control (Applied Biosystems)	
miRNA specific RT-qPCR assays	TaqMan MicroRNA assays (Applied Biosystems), miScript Primer Assays and miScript PCR Control Set (QIAGEN)	

## Results

Figure 1: Staining

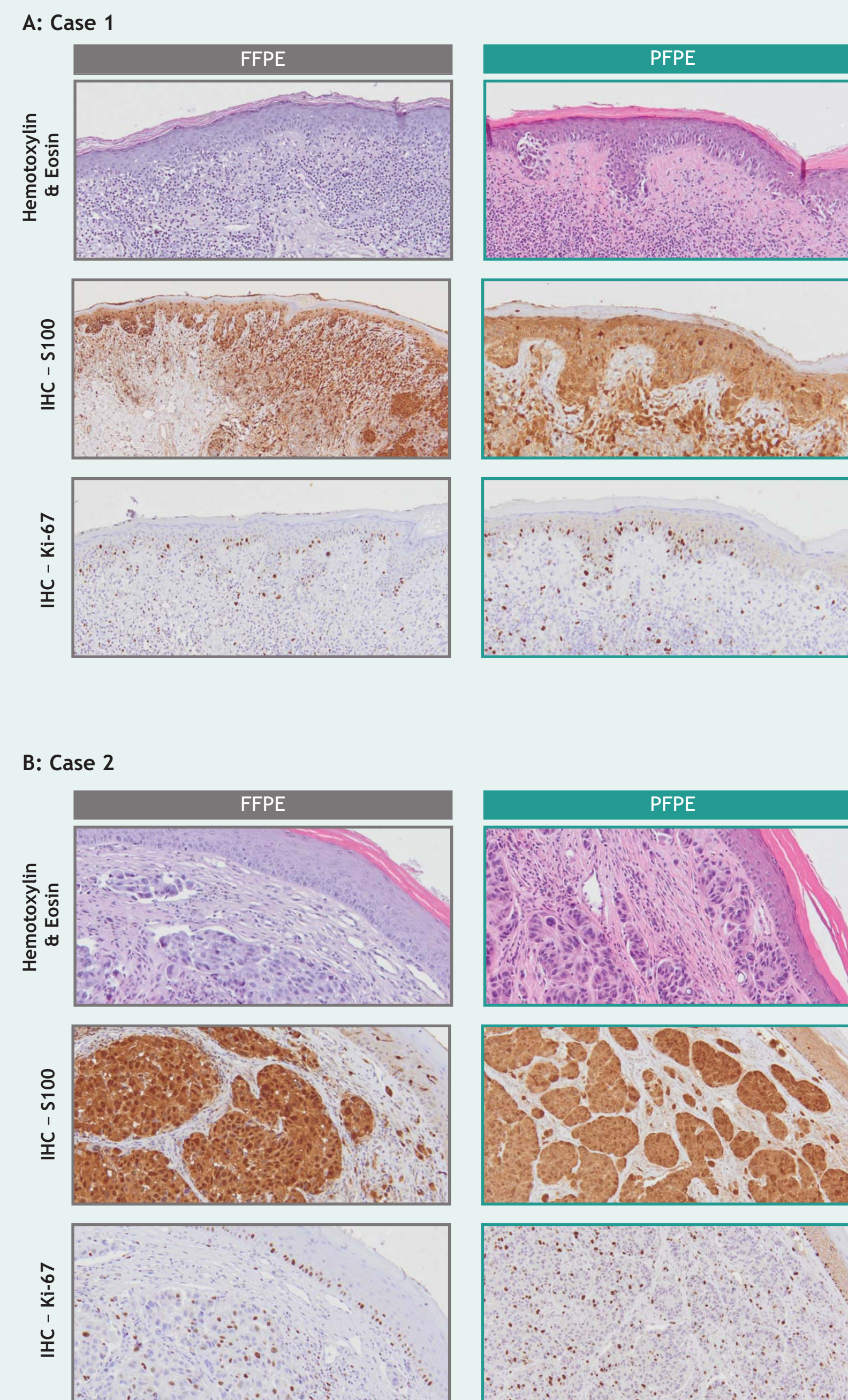


- Antibodies: Ki-67 clone MIB-1 (Dako M7240, 1:75 dilution); S100 polyclonal rabbit (Dako Z0311, 1:400 dilution)
- Epitope retrieval: FFPE 20 min at 98 °C in citrate buffer pH6, PFPE 10 min at 70 °C (Ki-67) or at 98 °C (S100) in Tris/EDTA buffer pH9

Figure 2: RNA from Cases 1-3, Purity and Integrity




Figure 3: mRNA Profiling with RNA from Cases 2 and 3

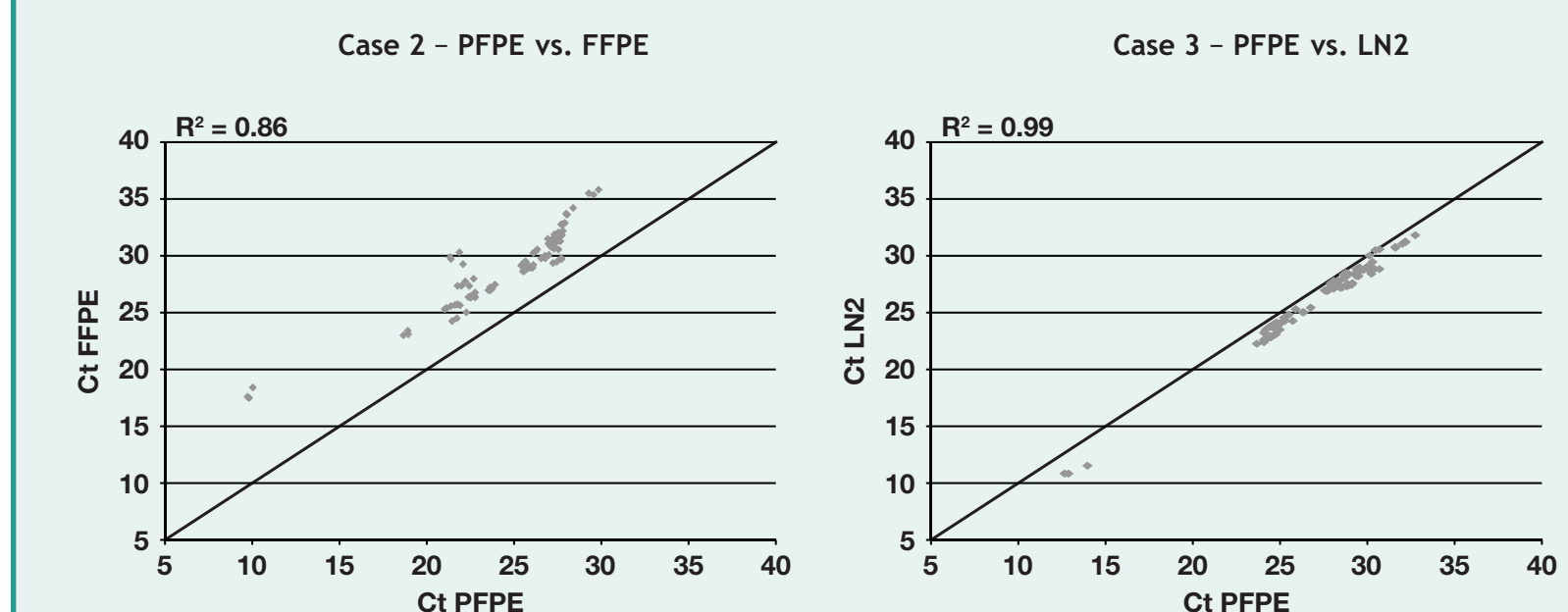

### Scatterplots for comparison of Ct values

- TaqMan Array: Human Endogenous Control with 32 control genes, plated in triplicates in a 96 well format
- Two Step RT-PCR with 2 µg RNA per plate
- R²: coefficient of determination

Figure 4: miRNA Profiling with miRNA from Cases 2 and 3

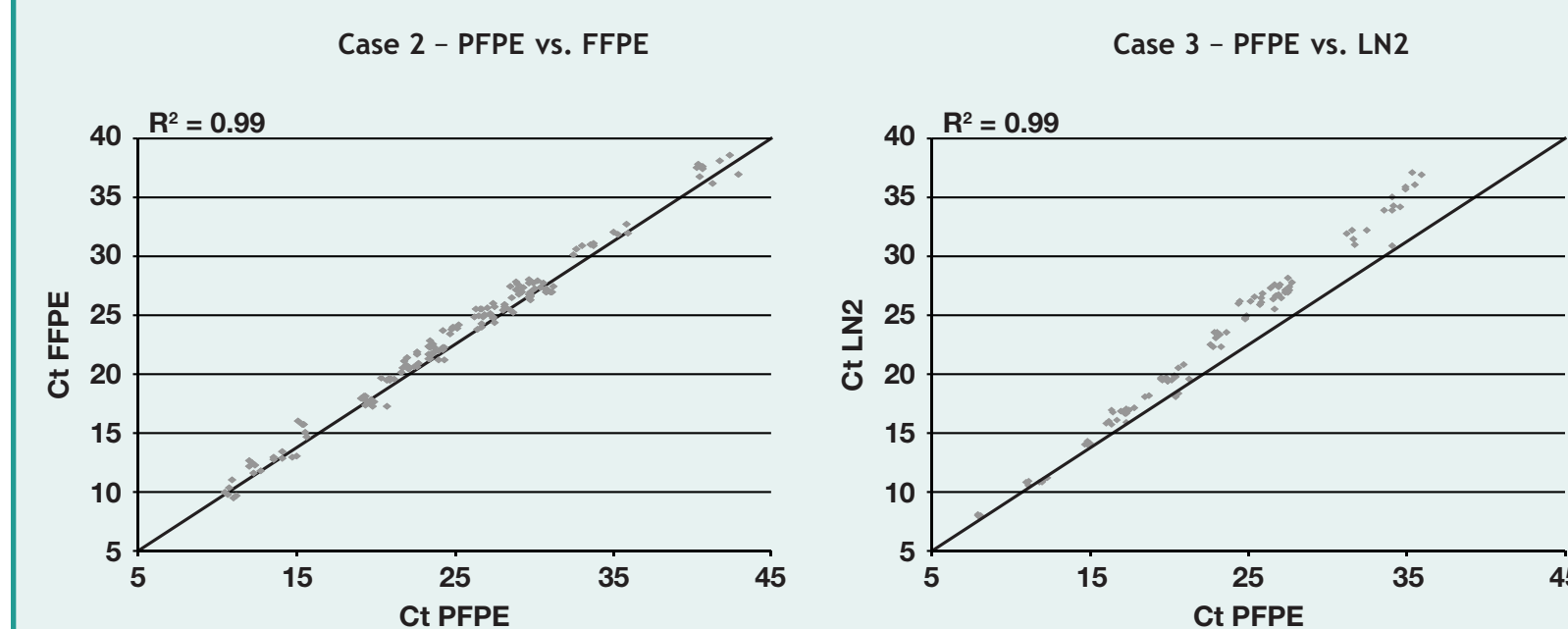

### Scatterplots with Ct values from 26 different noncoding small RNA specific RT-qPCR assays:

- 10x TaqMan MicroRNA Assays: miR10a, -16, -29a, -30b, -103, -137, -182, -192, let -7a, -7b
- 16x miScript miRNA Assays: miR9, -10a, -10b, -29a, -103, -125b, -143, -145, -155, -192
- 6x miScript PCR endogenous controls: RNUA1, RNU5A, RNU6B, SNORD25, SCARNA17, SNORA73A
- R²: coefficient of determination

## Summary

### Stabilization of human melanoma tissue with PAXgene Tissue:

- H&E Morphology: well preserved, comparable to FFPE
- Antigenicity: for Ki-67 and S100 comparable to FFPE with regard to location, staining intensity and number of stained cells
- Gene expression profile:
  - high concordance between RNA from PFPE and snap frozen tissue
  - low concordance between RNA from PFPE and FFPE
- Small, noncoding RNA profiles:
  - high concordance between RNA/miRNA from PFPE and snap frozen tissue
  - high concordance between RNA/miRNA from PFPE and FFPE

## Conclusions

This evaluation demonstrates the PAXgene Tissue System preserves morphology, antigenicity, RNA and miRNA profiles in human melanoma specimen

For Research Use Only. Not for use in diagnostic procedures. No claim or representation is intended to provide information for the diagnosis, prevention, or treatment of disease.