
April 2024

QIAxcel[®] Connect Safety Instructions and Quick-Start Guide

For use with QIAxcel ScreenGel[®] software v2.1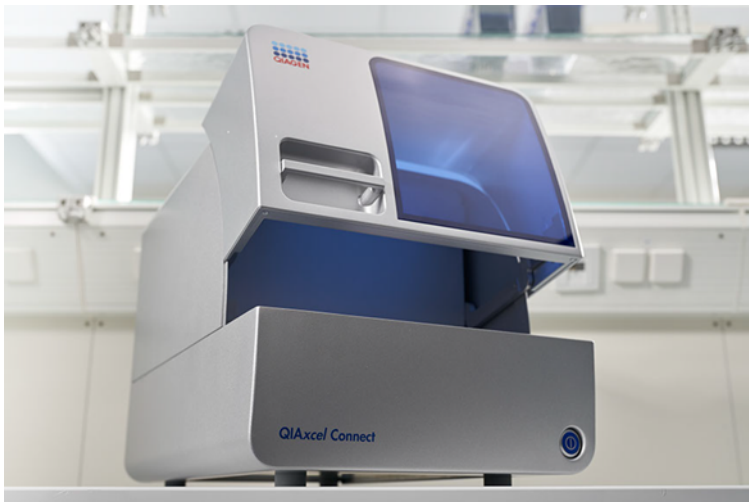

Sample to Insight

Table of Contents

Introduction	3
Safety Information	4
Proper use	6
Electrical safety	10
Environment	14
Chemicals	19
Waste disposal	21
Mechanical hazards	22
Installation and maintainance	23
Symbols on the QIAxcel Connect	27
Unpacking the QIAxcel Connect	30
Installation of the QIAxcel ScreenGel Operating Software	31
Installing the QIAxcel Connect	35
Using a QIAxcel Kit	38
Preparing the Buffer Tray	40
Preparing Markers	41
Loading the Buffer Tray	42
Installing the QIAxcel Gel Cartridge and Smart Key	44
Performing Cartridge Calibration	46
Running samples	48
Document Revision History	49

Introduction

This quick-start guide summarizes how to set up the QIAxcel Connect System. Before using the QIAxcel Connect System, it is essential that you read the QIAxcel Connect System User Manual carefully and pay particular attention to the safety information. The *QIAxcel Connect System User Manual* is available on the enclosed USB flash drive or under the **Resources** tab of the QIAxcel Connect product page (www.qiagen.com/qiaxcel-connect).

Safety Information

Before using the QlAxcel Connect, it is essential that you read this manual carefully and pay particular attention to the safety information. The instructions and safety information in the manual must be followed to ensure safe operation of the instrument and to maintain the instrument in a safe condition.

The following types of safety information appear throughout this manual:

WARNING

The term **WARNING** is used to inform you about situations that could result in personal injury to you or others.

Details about these circumstances are given in a box like this one.

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CAUTION

The term **CAUTION** is used to inform you about situations that could result in damage to the instrument or other equipment.

Details about these circumstances are given in a box like this one.

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The advice given in this manual is intended to supplement, not supersede, the normal safety requirements prevailing in the user's country.

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Improper use of the QIAxcel Connect may cause personal injuries or damage to the instrument.

The QIAxcel Connect must only be operated by qualified personnel who have been appropriately trained.

Servicing of the QIAxcel Connect must only be performed by QIAGEN® Field Service Specialists.

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Perform the maintenance as described in the Maintenance Procedures section of the *QIAxcel Connect System User Manual*. QIAGEN will charge for repairs required due to incorrect maintenance.

WARNING/ **Risk of personal injury and material damage**

CAUTION

The QIAxcel Connect is too heavy to be lifted by one person. To avoid personal injury or damage to the instrument, do not lift the instrument alone.

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WARNING **Risk of personal injury and material damage**

Do not attempt to move the QIAxcel Connect during operation.

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In case of emergency, switch off the QIAxcel Connect at the power switch at the front of the instrument and unplug the power cord from the power outlet.

CAUTION **Damage to the instrument**

If the Pressure 1 status is Low, increase system pressure before performing the unlatch command. Removing the cartridge and unlatching with reduced pressure may damage the instrument.

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CAUTION **Damage to the instrument**

Do not use bleach, solvents, or reagents containing acids, alkalis, or abrasives to clean the QIAxcel Connect.

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CAUTION **Damage to the instrument**

Avoid spilling water, chemicals, and liquids in general onto the QIAxcel Connect. Damage caused by water or chemical spillage will void your warranty.

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CAUTION **Damage to the instrument**

Do not stack instruments and do not place items on top of the QIAxcel Connect.

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Electrical safety

Disconnect the line power cord from the power outlet before servicing.

CAUTION Damage to electronics

Before switching on the instrument, make sure that the correct supply voltage is used. Use of incorrect supply voltage may damage the electronics. To check the recommended supply voltage, refer to the specifications indicated on the type plate of the instrument.

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WARNING

Electrical hazard

Use only the cable supplied with the QIAxcel Connect.

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WARNING

Electrical hazard

Any interruption of the protective conductor (earth/ground lead) inside or outside the instrument or disconnection of the protective conductor terminal is likely to make the instrument dangerous.

Intentional interruption is prohibited.

Lethal voltages inside the instrument

When the instrument is connected to line power, terminals may be live, and opening covers or removing parts is likely to expose live parts.

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To ensure satisfactory and safe operation of the QIAxcel Connect, follow the advice below:

- The line power cord must be connected to a line power outlet that has a protective conductor (earth/ground).
- Do not adjust or replace internal parts of the instrument.
- Do not operate the instrument with any covers or parts removed.
- If liquid has spilled inside the instrument, switch off the instrument, disconnect it from the power outlet, and contact QIAGEN Technical Services.
- When replacing the main fuse, replace only with the type and current rating specified on the type plate.

If the instrument becomes electrically unsafe, prevent other personnel from operating it, and contact QIAGEN Technical Services. The instrument may be electrically unsafe when:

- The line power cord appears to be damaged.
- It has been stored under unfavorable conditions for a prolonged period.
- It has been subjected to severe transport stresses.

Environment

Operating conditions

WARNING

Risk of explosion

The QIAxcel Connect is not designed for use in an explosive atmosphere.

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WARNING

Explosive atmosphere

The QIAxcel Connect is intended for use with reagents and substances supplied with the QIAGEN QIAxcel Kits. Use of other reagents and substances may lead to fire or explosion.

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CAUTION **Damage to the instrument**

Direct sunlight may bleach parts of the instrument and cause damage to plastic parts.

The QIAxcel Connect must be located out of direct sunlight.

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CAUTION Damage to the cartridge

When in use, the gel cartridge should not be removed from the Wash Park position of the buffer tray for more than 15 minutes. Failure to do so will cause the capillary tips to dry out, affecting correct function of the cartridge. Dry tips will void the warranty.

The capillary tips are made of glass and are very fragile. Take care not to knock the tips on any hard surfaces. Failure to do so will cause the capillary tips to break, affecting correct function of the cartridge. Broken tips will void the warranty.

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CAUTION **Damage to the cartridge**

If less than 12 samples are processed, fill the empty sample wells with the QX DNA Dilution Buffer or QX RNA Dilution Buffer. Failure to do so may cause damage to the capillary channels.

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CAUTION **Damage to the cartridge**

Direct sunlight may bleach the cartridge and reagents inside and cause damage to plastic parts.

The cartridge must be located out of direct sunlight.

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WARNING

Hazardous chemicals

Some chemicals used with this instrument may be hazardous or may become hazardous after completion of purification.

Always wear safety glasses, gloves, and a lab coat.

The responsible body (for example, a laboratory manager) must take the necessary precautions to ensure that the surrounding workplace is safe, and that the instrument operators are suitably trained and not exposed to hazardous levels of infectious agents as defined in the applicable Material Safety Data Sheets (MSDSs) or the OSHA^{1*}, ACGIH[†], or COSHH[‡] documents.

Venting for fumes and disposal of waste must be in accordance with all national, state, and local health and safety regulations and laws.

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* OSHA – Occupational Safety and Health Organization (United States of America)

† ACGIH – American Conference of Government Industrial Hygienists (United States of America)

‡ COSHH – Control of Substances Hazardous to Health (United Kingdom)

WARNING**Risk of fire**

When cleaning the QIAxcel Connect with alcohol-based disinfectant, leave the QIAxcel Connect doors open to allow flammable vapors to disperse.

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Waste disposal

Used labware and containers may contain hazardous chemicals. Such wastes must be collected and disposed of properly according to local safety regulations.

For information on how to dispose of the QIAxcel Connect, see Appendix A of the *QIAxcel Connect System User Manual*.

Mechanical hazards

The cartridge door and sample door of the QIAxcel Connect must remain closed during operation of the instrument.

WARNING

Moving parts

To avoid contact with moving parts during operation of the QIAxcel Connect, the instrument must be operated with the cartridge door and sample door closed.

If the door sensors are not functioning correctly, contact QIAGEN Technical Services.

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WARNING Risk of overheating

To ensure proper ventilation, maintain a minimum clearance of 10 cm (3.94 in.) at the sides and rear of the QIAxcel Connect.

Slits and openings that ensure the ventilation of the QIAxcel Connect must not be covered.

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**WARNING/
CAUTION**

Risk of personal injury and material damage

The covers of the QIAxcel Connect must only be removed by qualified personnel appropriately trained by QIAGEN.

Servicing of the QIAxcel Connect must only be performed by QIAGEN-authorized representatives.

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WARNING

Risk of fire

Do not allow cleaning fluid or decontamination agents to come into contact with the electrical parts of the QIAxcel Connect. Take special care while cleaning the sample plate drawer that no liquid is spilled into the inside of the instrument.

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WARNING

Risk of electric shock

Do not open any panels on the QIAxcel Connect other than described in this user manual.

Risk of personal injury and material damage

Only perform maintenance that is specifically described in this user manual.

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WARNING**Risk of explosion**

When cleaning the QlAxcel Connect with alcohol-based disinfectant, allow flammable vapors to disperse.

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WARNING**Risk of fire or explosion**

When using ethanol or ethanol-based liquids on the QlAxcel Connect, handle such liquids carefully and in accordance with the required safety regulations. If liquid has been spilled, wipe it off and allow flammable vapors to disperse.

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Unpacking the QIAxcel Connect

1. Identify a suitable place for the QIAxcel Connect on the laboratory bench (see Table 1 below).
2. Ensure that the bench is dry, clean, and vibration-proof, and has additional space for accessories.
3. Unpack and place the QIAxcel Connect onto the bench.

Table 1. QIAxcel Connect specifications and operating conditions

QIAxcel Connect specifications

Dimensions (door closed)	Width: 372 mm (14.64 in.) Height: 408 mm (16.06 in.) Depth: 572 mm (22.44 in.)
Dimensions (doors open)	Width: 372 mm (14.64 in.) Height: 599 mm (23.58 in.) Depth: 572 mm (22.44 in.)
Mass	32 kg (70.55 lb.)
Operating conditions	
Air temperature	15–30°C (59–86°F)
Relative humidity	10–75% (noncondensing)
Altitude	Up to 2000 m (6500 ft.)
Power	100–240 V AC, 50–60 Hz, 100 VA

Installation of the QIAxcel ScreenGel Operating Software

The QIAxcel ScreenGel software v2.1 or higher must be installed before connecting the QIAxcel Connect to the computer for the first time. The QIAxcel ScreenGel software can be installed from the USB drive supplied with the QIAxcel Connect. The QIAxcel ScreenGel software version lower than 2.0 will not work with the QIAxcel Connect.

Things to do before starting

1. If you are using a computer other than provided with the QIAxcel Connect, ensure that the computer meets the minimum requirements (see *QIAxcel Connect System User Manual*).
2. Use a Windows® account with administrator privileges and close any programs running on the computer.
3. Make sure that the QIAxcel Connect is not connected to the computer (i.e., unplug the cable from the computer).
4. The installation of the QIAxcel ScreenGel software is possible via the supplied USB drive. Internet access is required for the software download from the product page at www.qiagen.com/qiaxcel-connect.

Note: We recommend having no other software installed on the computer for operating the QIAxcel Connect other than the QIAxcel ScreenGel software.

5. Insert the enclosed USB drive in the computer's USB port.
6. Double-click **My Computer**, and double-click the **USB drive**. Launch the **QIAxcelSetup.exe** program. This will start the QIAxcel ScreenGel software installation. The setup wizard will guide you through the setup process.

Note: It is possible to switch the language of the installed software later using the File menu.

Software installation

1. To start the actual installation procedure, select the **Install QIAXcel ScreenGel Software** option.

Note: To read the user manual in the selected language, select the **User Manual** option.

Note: The **Microsoft Visual C++ 2015-2022 x86 Redistributable** will be installed automatically during QIAXcel ScreenGel wizard installation.

2. ScreenGel software v2.1 requires version 4.7.2 of the Microsoft .NET Framework. This .NET version is already installed in Windows 10 version 1803 or higher. To use ScreenGel software v2.1 with an earlier version of Windows 10, the Microsoft® .NET Framework 4.7.2 must be installed manually.
3. The QIAXcel ScreenGel setup wizard opens. Click **Next**.
4. Accept the license agreement and click **Next**.
5. Select the program installation path. The default path is `c:\Program Files (x86)\QIAGEN\QIAXcel ScreenGel 2\`. Click **Next**.

Note: If `c:\Program Files (x86)\QIAGEN\QIAXcel ScreenGel 2\` is selected on a Windows 64 bit system, the installation will appear in `c:\Program Files\x86\QIAGEN\QIAXcel\ScreenGel`.

6. Select the data path for acquired data and all other application data. The default path is `c:\ProgramData\QIAGEN\QIAXcel\ScreenGel 2\`. Click **Next**.

Note: It is not possible to choose `c:\Program Files (x86)\QIAGEN\QIAXcel ScreenGel 2\` also as the data path. It is also not possible to choose the system root path (e.g., `c:\windows`) as the data path.

Note: All paths in this manual refer to the data path as `%DATA_DIR%`.

Note: The `%DATA_DIR%` directory and its subdirectories can be opened from the Windows Explorer directly from QIAXcel ScreenGel using the **File/Open Data Directory** menu item.

7. To begin the installation, click **Install**. This may take some time.
8. The QIAGEN QIAxcel Connect Driver Installer dialog opens, if the driver is not yet installed on the computer. To install the driver, click **Next**. To complete the driver installation, click **Finish**.
9. To complete the QIAxcel ScreenGel setup wizard, click **Finish**.
10. To close the installation window, click **Exit**.
11. After the installation, the QIAxcel ScreenGel can be launched either from the Windows Start menu under **QIAGEN/QIAxcel** or from the desktop icon.

First log in

1. For first log in, log in as **Administrator** and click **Login**.

Note: For this first login, no password is required.

QIAxcel ScreenGel 2.1.0.12

QIAGEN

Mode
DNA

User ID
Administrator

Password

Cancel

Login

Sample to Insight

Figure 1. First Login.

2. Provide a valid password for the Administrator account. Leave the field for the old password empty, then provide a new valid password and confirm it by clicking **OK** (see Figure 2).

Note: The password must contain one uppercase character, one lowercase character, and one digit. The minimum length of the password is eight characters.

QIAGEN Password for user Administrator has expired.
Enter a new password.

Re-enter your old password

Enter a new password

Confirm the new password

OK Cancel

Figure 2. Creating the password for the Administrator.

3. The Configuration environment is displayed by default. Select the **User Manager** tab.
4. Create user accounts for all relevant users. For more information, refer to the User Management section of the *QIAXcel Connect System User Manual*.
5. Configure which COM port is to be used for the connection to the QIAXcel Connect system (default adjustment is COM1) and the global settings for the QIAXcel ScreenGel software. For more detailed information, refer to the Settings section of the *QIAXcel Connect System User Manual*.
6. Configure the global settings for the QIAXcel ScreenGel software. For further details, refer to the Settings section of the *QIAXcel Connect System User Manual*.

Installing the QIAxcel Connect

The following steps must be performed before the QIAxcel Connect can be operated:

1. Make sure that the QIAxcel Connect has equilibrated to ambient temperature before operating it and that the power switch of the QIAxcel Connect is in the **Off** position.
2. Open the sample door.
3. Remove the transport lock by turning counterclockwise (Figure 3 on the next page). Save the lock for future use. A storage bay for the transport lock is located under the Service Door for storage use (at the top of the instrument).
4. Install the N₂ cylinder (Figure 4 on the next page). Screw the N₂ cylinder clockwise into the N₂ cylinder port (to the right of the sample plate holder). Turn until the needle inside the port pierces the N₂ cylinder. The cylinder should only be finger tight; do not overtighten. Gently push the N₂ cylinder down until it is in the stowed (down) position.

Important: Use only N₂ cylinders provided by QIAGEN (cat. no. 929705). Alternatively, external, clean, non-condensing compressed nitrogen (min. 50 psi, max. 75 psi) can be connected to the external N₂ port (Figure 5) via the supplied urethane tubing.

5. Connect the enclosed USB cable to the USB port of the QIAxcel Connect (Figure 5) and the computer.
6. Connect the power cord to the QIAxcel Connect power connection (Figure 5) and into a grounded power outlet.

Figure 3. Removing the transport lock.

Figure 4. Buffer tray holder (a), sample plate holder (b), and N2 port (c) location.

Figure 5. Power connection (a), external N2 (b), and USB port (c).

Using a QIAxcel Kit

1. Ensure all equipment and reagents are equilibrated to room temperature (15–25°C) before use.
2. Add 10 mL QX (HS) Wash Buffer to both reservoirs of the QX Cartridge Stand (provided with the instrument) and cover with 2 mL mineral oil.
3. Remove the QIAxcel Gel Cartridge from its packaging and carefully wipe off any soft gel debris from the capillary tips using a soft tissue.

Note: The QIAxcel Gel Cartridge is provided together with a smart key, which is used for fully automatic cartridge detection. Do not remove the smart key from the cartridge.

4. Remove the purge cap seal from the back of the QIAxcel Gel Cartridge (Figure 6) and place it in the QX Cartridge Stand (Figure 7). Retain the purge port seal in case you need to store the QIAxcel Gel Cartridge.

Note: Use a soft tissue to wipe off any gel that may have leaked from the purge port.

Note: Ensure that the capillary tips are submerged in QX Wash Buffer.

5. **Important:** Cartridges must be allowed to equilibrate at room temperature for at least 20 min prior to use. Place the gel cartridge in the QX Cartridge Stand, protected from light with the QX Cartridge Stand Cover.

Note: Incorrect cartridge handling or shorter equilibration times may cause damage to your QIAxcel Connect instrument and could result in a loss of warranty.

Figure 6. Preparing the QIAxcel Gel Cartridge.

Figure 7. Incubating in the QX Cartridge Stand.

Preparing the Buffer Tray

1. Allow all reagents to equilibrate to room temperature (15–25°C) before use.
2. Wash the buffer tray (Figure 8) with warm water and rinse thoroughly with deionized water.
3. Fill the WP (Wash Park) and WI (Wash Inject) positions of the buffer tray with 8 mL QX (HS) Wash Buffer.
4. Fill the BUFFER position of the buffer tray with 18 mL QX Separation Buffer, QX HS Separation Buffer, or QX FA Separation Buffer, depending on the kit in use.
5. Carefully add mineral oil to cover all 3 positions to prevent evaporation: 2 mL mineral oil to the WP and WI positions and 4 mL mineral oil to the BUFFER position.

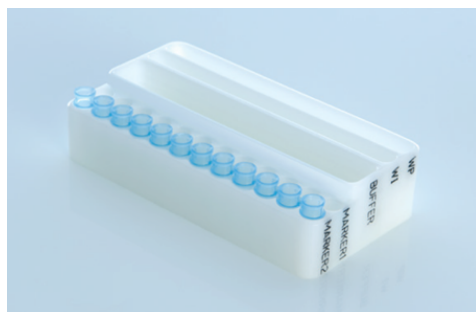
Figure 8. Preparing the buffer tray.

Preparing Markers

1. Load 15 μL QX (HS) Alignment Marker into each tube of a QX 0.2 mL 12-Tube Strip. Add 1 drop of mineral oil to each tube and place the strip into the MARKER1 position of the buffer tray.
2. Load 15 μL QX (HS) Intensity Calibration Marker into each tube of a QX Colored 0.2 mL 12-Tube Strip. Make sure there are no air bubbles, add a drop of mineral oil to each tube, and insert the strip into the MARKER2 position of the buffer tray.

Loading the Buffer Tray

1. Close the cartridge door and the sample door.

Note: The QIAxcel Connect's cartridge and sample doors must remain closed during operation of the instrument. Opening either door during operation will cause the system to stop any action it is currently performing.

2. Switch on the QIAxcel Connect at the power switch. Switch on the computer and launch the QIAxcel ScreenGel software from the Windows Start menu under **QIAGEN/QIAxcel ScreenGel** or from the desktop icon.
3. Select a mode (DNA or RNA) and log in.

Note: The  icon indicates that the connection is being established, and the  icon shows that the QIAxcel Connect is connected. In the case that the instrument could not be connected, a message will be shown notifying that the instrument is not available. If you did not switch on the instrument at the moment, click **Instrument not needed**. If you need the instrument, click **Troubleshoot**. Follow the instructions in the message. For more detailed instructions, refer to the System setup subsection under the Troubleshooting section, within the *QIAxcel Connect System User Manual*. Close the message. To retry to connect to the instrument, click the  icon.

4. Click  in the Status Information panel at the left to move the buffer tray holder to the front of the instrument. Wait until the buffer tray holder has reached its stop position.
5. Open the sample door and carefully place the filled buffer tray into the buffer tray holder, making sure to place the buffer tray in the correct orientation (Figure 4).

Note: Be careful not to spill any solutions in the instrument or cause any cross-contamination between buffers loaded on the buffer tray.

Note: The 12-tube strips should be positioned towards the front of the instrument with the buffers towards the back.

Important: Always remove the buffer tray before moving the QIAxcel Connect.

6. Close the cartridge and the sample door and click the  icon in the Status Information panel to move the buffer tray holder to the Wash Park position.

Note: Leave the sample door open if you want to load your samples immediately afterward.

Note: If you close the sample door, the buffer tray automatically moves to the Wash Park position after a period of 5 minutes.

Installing the QIAxcel Gel Cartridge and Smart Key

1. Remove the QIAxcel Gel Cartridge from the QX Cartridge Stand.
2. Open the cartridge door and place the QIAxcel Gel Cartridge into the QIAxcel Connect. The cartridge description label should face the front and the purge port should face the back of the instrument (Figure 9).

Note: Ensure the purge cap seal has been removed as described above.

3. Insert the smart key into the smart key socket near the cartridge (Figure 10). The smart key can be inserted in either direction.

Note: The system will not recognize the cartridge and will not operate if the smart key is not inserted.

4. Close the cartridge door.

Figure 9. Installation of QIAxcel Gel Cartridge.

Figure 10. Insertion of smart key.

Performing Cartridge Calibration

The QIAxcel Gel Cartridges must be calibrated before they are used for the first time. Every new cartridge requires an intensity calibration before the first sample analysis. The calibration normalizes the intensities of each capillary by applying a factor for every subsequent run with this cartridge. See the *Calibrating a cartridge* section in the *QIAxcel Connect System User Manual* for more details.

1. Open the **Calibration** tab of the Service environment. Here, click **Start Calibration** to launch a calibration run.

Note: The total running time for the calibration process is approximately 16 minutes.

2. In the next dialog, confirm that the QX Intensity Calibration Marker is loaded. Optionally, enter the lot ID (refer to Figure 11). Start the calibration by clicking **OK**.

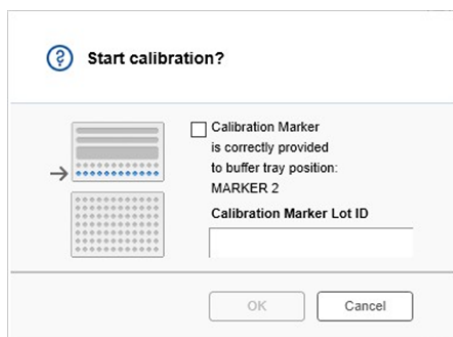
Figure 11. Start calibration dialog window.

3. Once the calibration process is complete, the calibration results are displayed next to the gel image or in the electropherogram view (Figure 12). The result table shows the area, the calibration factor, and the result (Pass or Fail) for each capillary channel.

4. When the calibration process is completed successfully, click **Finish Calibration** to accept the calibration result. The cartridge is now calibrated.

Note: If one or more capillary channels fail, repeat the calibration process and overwrite the previous results. If the issue persists, contact QIAGEN Technical Services.

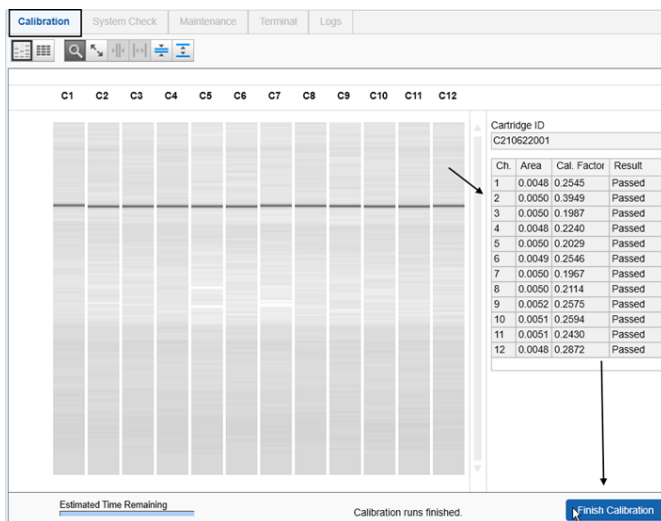
Figure 12. Calibration results of a QIAxcel Gel Cartridge.

Running samples

Refer to the Running a process section in the *QIAxcel Connect System User Manual* for detailed instructions.

Document Revision History

Date	Changes
10/2021	Initial version of the QIAxcel Connect Quick Start Guide
04/2024	Update the QIAxcel Connect Quick Start Guide according to version 2.1 of the QIAxcel Connect ScreenGel software.

For further support with technical difficulties, please contact QIAGEN Technical Services at www.qiagen.com

For up-to-date licensing information and product-specific disclaimers, see the respective QIAGEN user manual. QIAGEN kit handbooks and user manuals are available at www.qiagen.com or can be requested from QIAGEN Technical Services or your local distributor.

Trademarks: QIAGEN®, Sample to Insight®, QIAxcel®, ScreenGel® (QIAGEN Group); Microsoft®, Windows® (Microsoft Corporation). Registered names, trademarks, etc. used in this document, even when not specifically marked as such, are not to be considered unprotected by law.

HB-294-002 04/2024 © 2024 QIAGEN, all rights reserved.

Page left intentionally blank.