

DNeasy[®] Plant Mini Kit

The DNeasy Plant Mini Kit (cat. nos. 69104 and 69106) can be stored at room temperature (15–25°C) for up to 1 year if not otherwise stated on label.

Further information

- *DNeasy Plant Handbook*: www.qiagen.com/HB-1166
- Safety Data Sheets: www.qiagen.com/safety
- Technical assistance: support.qiagen.com

Notes before starting

- Perform all centrifugation steps at room temperature (15–25°C).
- If necessary, redissolve any precipitates in Buffer AP1 and Buffer AW1 concentrates.
- Add ethanol to Buffer AW1 and Buffer AW2 concentrates.
- Preheat a water bath or heating block to 65°C.

1. Disrupt samples (≤ 100 mg wet weight or ≤ 20 mg lyophilized tissue) using the TissueRuptor[®], the TissueLyser II or a mortar and pestle.
2. Add 400 μ l Buffer AP1 and 4 μ l RNase A. Vortex and incubate for 10 min at 65°C. Invert the tube 2–3 times during incubation.

Note: Do not mix Buffer AP1 and RNase A before use.

3. Add 130 μ l Buffer P3. Mix and incubate for 5 min on ice.
4. **Recommended:** Centrifuge the lysate for 5 min at 20,000 x g (14,000 rpm).
5. Pipet the lysate into a QIAshredder spin column placed in a 2 ml collection tube. Centrifuge for 2 min at 20,000 x g.

6. Transfer the flow-through into a new tube without disturbing the pellet if present. Add 1.5 volumes of Buffer AW1, and mix by pipetting.
7. Transfer 650 μ l of the mixture into a DNeasy Mini spin column placed in a 2 ml collection tube. Centrifuge for 1 min at $\geq 6000 \times g$ (≥ 8000 rpm). Discard the flow-through. Repeat this step with the remaining sample.
8. Place the spin column into a new 2 ml collection tube. Add 500 μ l Buffer AW2, and centrifuge for 1 min at $\geq 6000 \times g$. Discard the flow-through.
9. Add another 500 μ l Buffer AW2. Centrifuge for 2 min at $20,000 \times g$.
Note: Remove the spin column from the collection tube carefully so that the column does not come into contact with the flow-through.
10. Transfer the spin column to a new 1.5 ml or 2 ml microcentrifuge tube.
11. Add 100 μ l Buffer AE for elution. Incubate for 5 min at room temperature (15–25°C). Centrifuge for 1 min at $\geq 6000 \times g$.
12. Repeat step 11.



Scan QR code for handbook.

For up-to-date licensing information and product-specific disclaimers, see the respective QIAGEN kit handbook or user manual.

Trademarks: QIAGEN[®], Sample to Insight[®], DNeasy[®], TissueRuptor[®] (QIAGEN Group). 1101205 03/2016 HB-0542-003 © 2016 QIAGEN, all rights reserved.